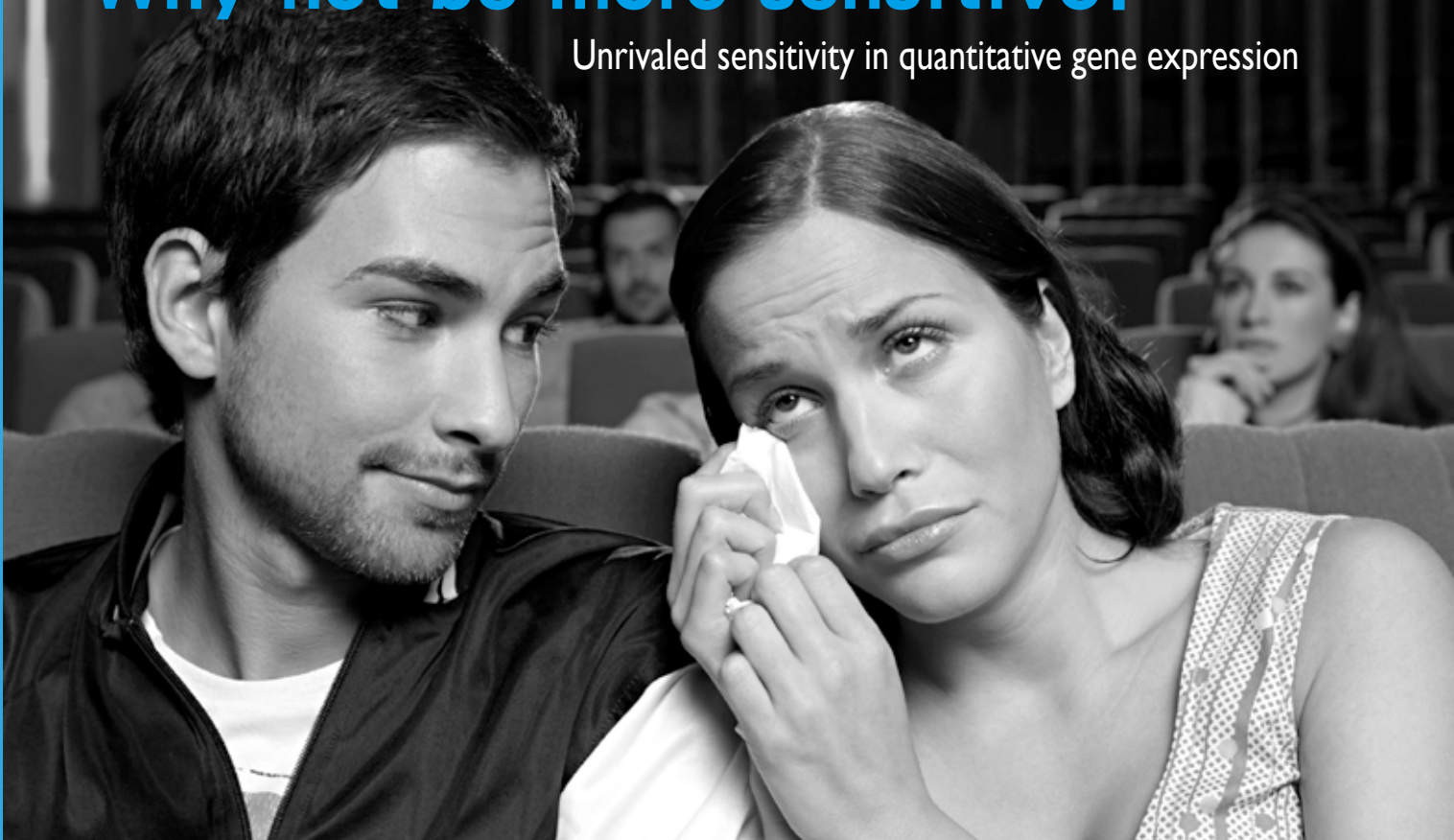
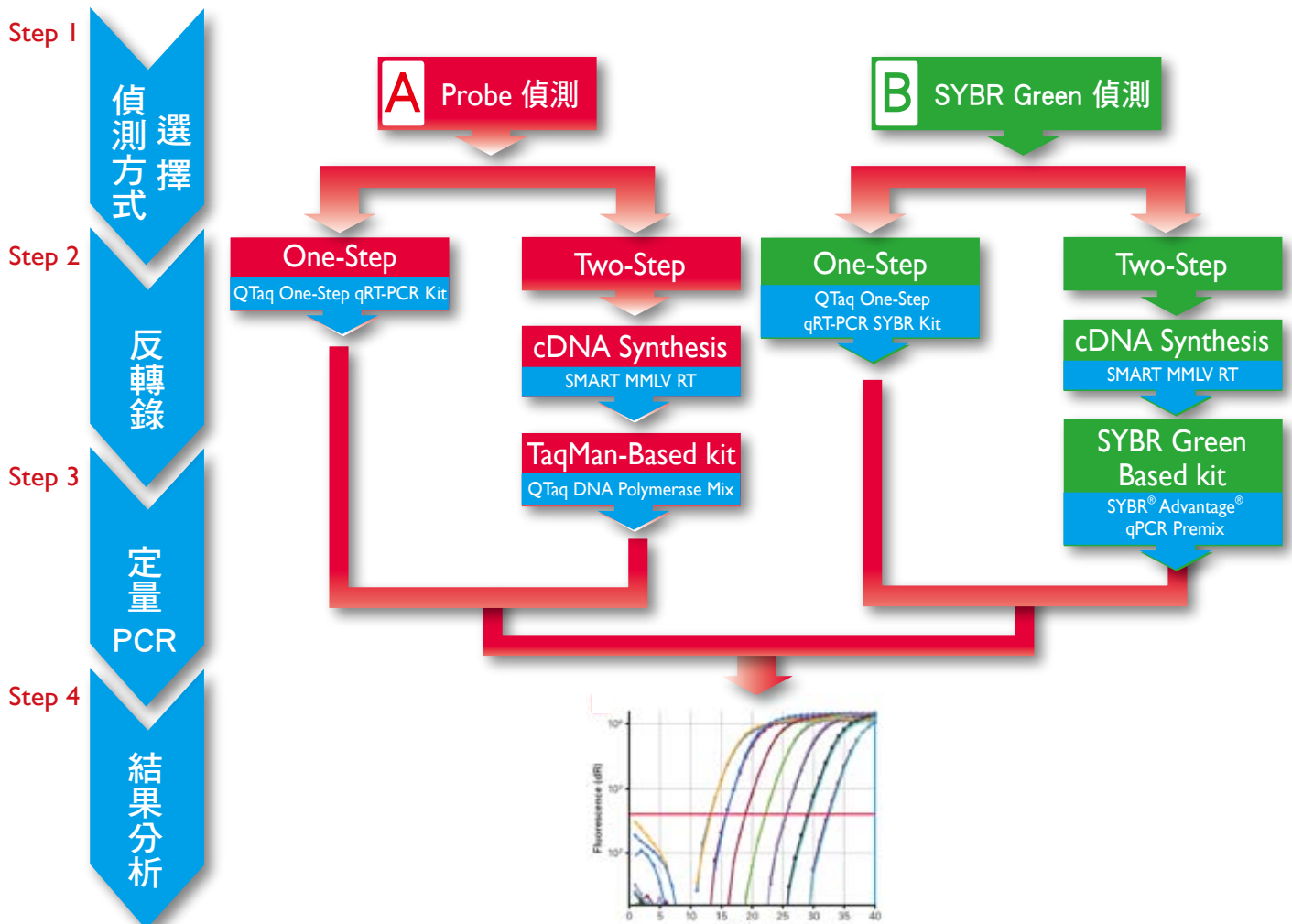
Why not be more sensitive?

Unrivalled sensitivity in quantitative gene expression

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• ROUND 1 • 專一性 (Specificity) 測試:

START!!

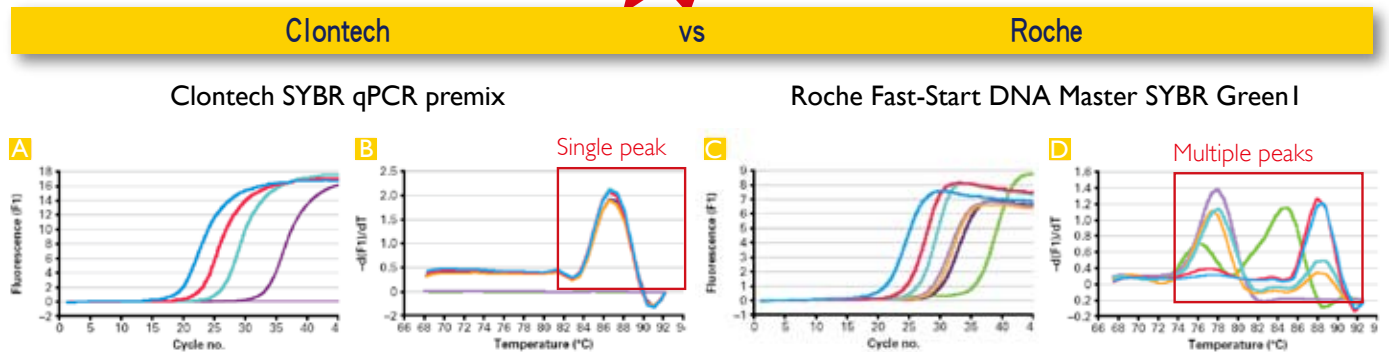
Figure 1. Reaction specificity—performance of the Clontech SYBR Advantage qPCR Premix vs. the Roche Fast Start DNA Master SYBR Green I using a Roche LightCycler. The results for the Clontech reagent are shown in Panels A and B, and those for the Roche reagent are shown in Panels C and D. The cycling conditions for the Clontech reagent consisted of 1 cycle at 95°C for 10 sec, followed by 45 cycles at 95°C for 5 sec and 60°C for 20 sec. For the Roche reagent, the cycling conditions consisted of 1 cycle at 95°C for 10 min, followed by 45 cycles at 94°C for 10 sec, 55°C for 5 sec, and 72°C for 10 sec.

• ROUND 2 • 專一性 (Specificity) 測試:

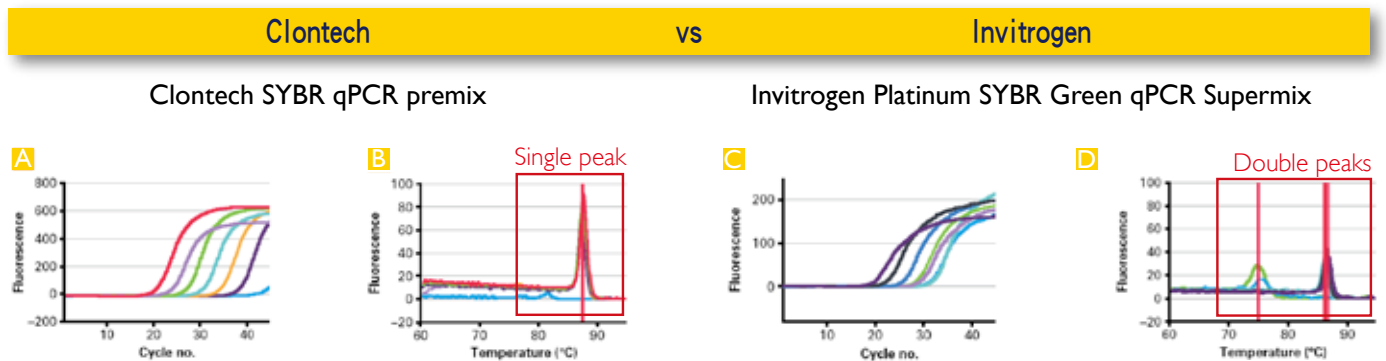
Figure 2. Reaction specificity—performance of the Clontech SYBR Advantage qPCR Premix vs. the Invitrogen Platinum SYBR Green qPCR Supermix UDG using a Cepheid Smart Cycler. The results for the Clontech reagent are shown in Panels A and B, and those for the Invitrogen reagent are shown in Panels C and D. The cycling conditions for the Clontech reagent consisted of 1 cycle at 95°C for 2 min, followed by 45 cycles at 95°C for 5 sec and 60°C for 20 sec. For the Invitrogen reagent, the cycling conditions consisted of 1 cycle at 95°C for 2 min, followed by 45 cycles at 95°C for 15 sec and 60°C for 30 sec.

• ROUND 3 • 效能(Efficiency)測試

Clontech

vs

ABI

Clontech SYBR qPCR premix

ABI SYBR Green PCR Master Mix

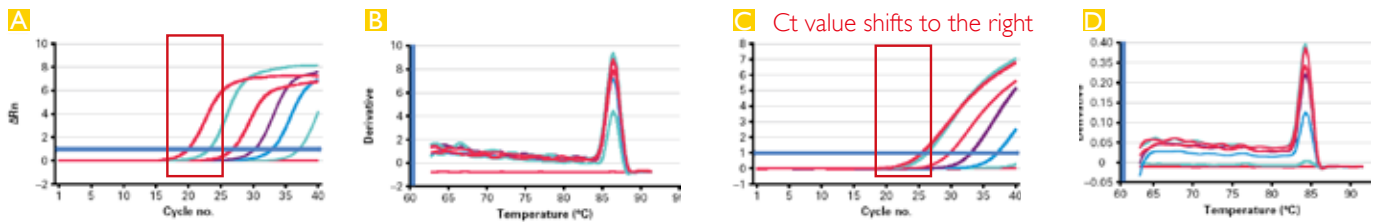
Figure 3. Amplification efficiency—performance of the Clontech SYBR Advantage qPCR Premix vs. the Applied Biosystems SYBR Green PCR Master Mix using an ABI PRISM 7000 Sequence Detection System. The results for the Clontech reagent are shown in Panels A and B, and those for the Applied Biosystems reagent are shown in Panels C and D. The cycling conditions for the Clontech reagent consisted of 1 cycle at 95°C for 10 sec, followed by 40 cycles at 95°C for 5 sec and 60°C for 31 sec. For the Applied Biosystems reagent, the cycling conditions consisted of 1 cycle at 95°C for 10 min, followed by 45 cycles at 95°C for 15 sec and 60°C for 1 min.

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Clontech April 2006 14-15

www.clontech.com

SYBR Advantage qPCR Premix Outperforms Competitors

Clearly, the Clontech SYBR Advantage qPCR Premix outperforms much of the competition in terms of specificity (single product as judged by melting curves), efficiency (amplification plots [shown] and standard curves [not shown]) and sensitivity (lower Ct values). Moreover, the overall

time for the qPCR analysis can be shortened dramatically—nearly in half compared to the times for other qPCR chemistries—when using the Clontech SYBR Advantage qPCR Premix.

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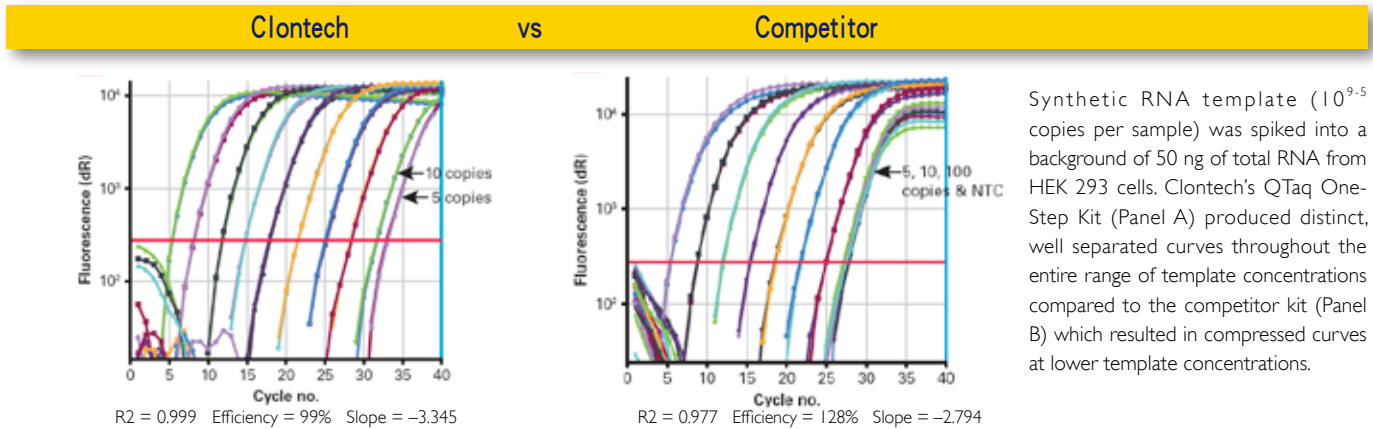
Synthetic RNA template (10^{9-5} copies per sample) was spiked into a background of 50 ng of total RNA from HEK 293 cells. Clontech's QTag One-Step Kit (Panel A) produced distinct, well separated curves throughout the entire range of template concentrations compared to the competitor kit (Panel B) which resulted in compressed curves at lower template concentrations.

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
Denatured time (min) 1 5

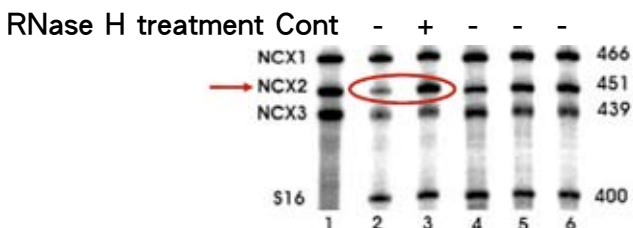Denatured temperature 94 97 100

"We show that the template of one gene of a multigene family becomes more readily available for PCR amplification only after RNase H treatment, while two other genes are not affected by RNase H treatment."

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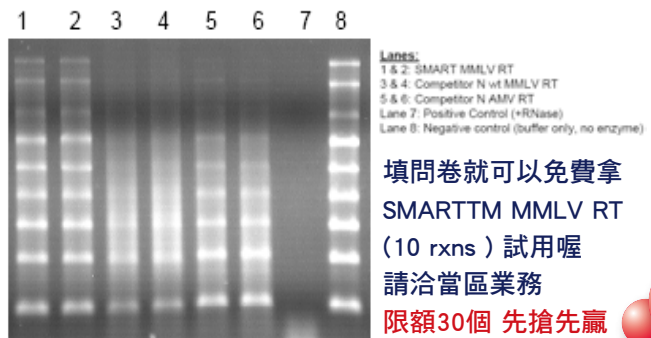
Figure 1. SMART MMLV RT has less RNase activity than the competition. An RNA Ladder was incubated with 200 units of either SMART MMLV RT (Lanes 1 & 2), or another supplier's MMLV RT (Lanes 3 & 4) and AMV RT (Lanes 5 & 6) at 37°C for 1 hr; and then analyzed by formaldehyde/agarose gel electrophoresis. Lane 7: positive control; Lane 8: negative control (buffer only, no enzyme).

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