Mir-X™ MicroRNA Quantification

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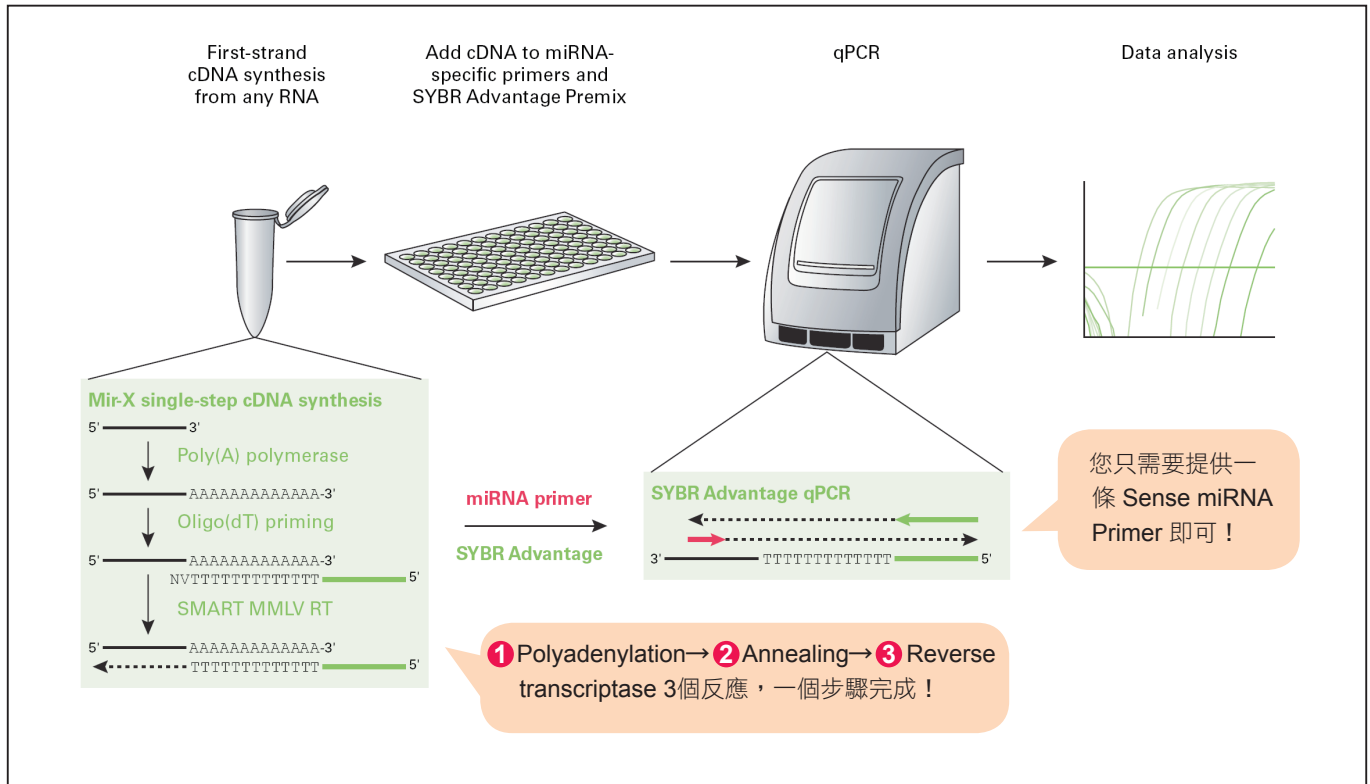
<http://microrna.sanger.ac.uk/sequences/>

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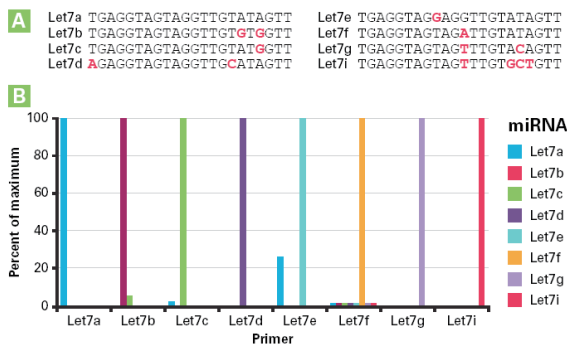
Figure 1. Specific quantification of Let7 miRNA variants. Using miRNA-specific primers (Panel A), Mir-X qRT-PCR was able to specifically detect and quantify each member of a series of 8 synthetic Let7 variants that had been spiked into a background of yeast polyA+ RNA (Panel B). The primers detected each of their corresponding Let7 miRNA cognates, but did not detect the off-target variants in 63 of 64 possible combinations.

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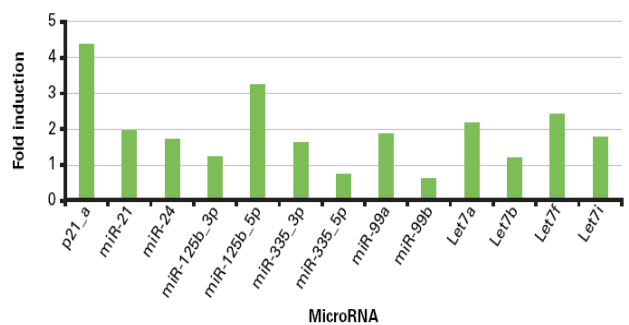
Figure 2. Trichostatin A treatment alters miRNA expression in mouse ES cells. Mouse embryonic stem cells were harvested either prior to or after being treated with trichostatin A (TSA) for 18 hr. RNA was prepared from the cells, and was then analyzed by Mir-X miRNA qRT-PCR using primers specific for the 12 indicated miRNAs and for a p21 control mRNA known to be induced by TSA.

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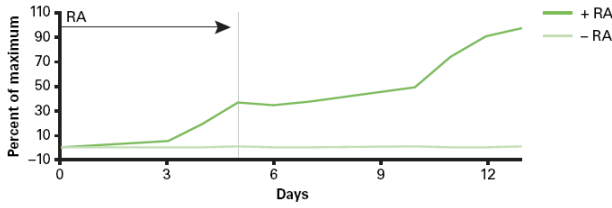
Figure 3. Induction of miR-9 in mouse P19 cells.

P19 mouse embryonal carcinoma cells were plated on agarose-coated petri dishes and allowed to form embryoid bodies (EB) in the presence or absence of retinoic acid (RA). After five days of culture, EBs were dissociated with trypsin and replated on tissue culture dishes without RA. Cells were harvested on days 3–7 and 10–13, and the induction of miR-9 was followed and quantified using the Mir-X miRNA qRT-PCR SYBR Kit and primers specific for miR-9 and U6 (as a normalization control).

Kimberly-Clark PROFESSIONAL **Kimcare**

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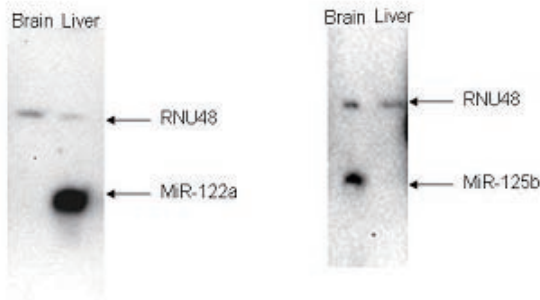
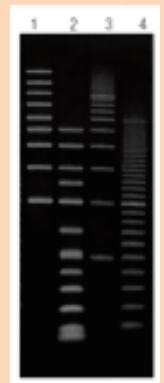
Ordering information

貨號	品名	包裝
NB-0001	miRNA northern blot kit with gel	For 2 miRNA probes each for monitoring 14 RNA samples or 6 miRNA probes each for 2 samples
NB-0002	miRNA northern blot kit without gel	For 2 miRNA probes each for monitoring 14 RNA samples or 6 miRNA probes each for 2 samples
MP-0XXX	miRNA probe	20 rxn