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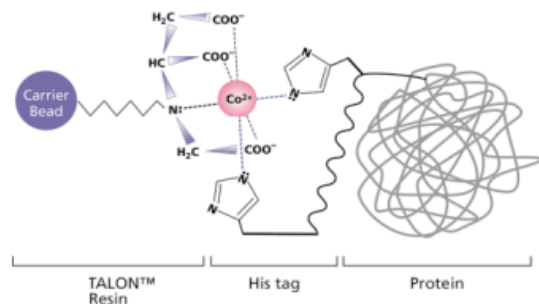
- Crystallography
- Functional Assay
- Structural investigation

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

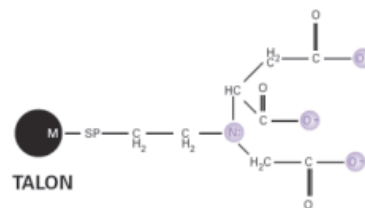
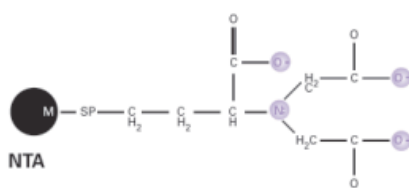
Two common chelating ligands used in MACS

TALON 能專一結合 His-tagged 蛋白

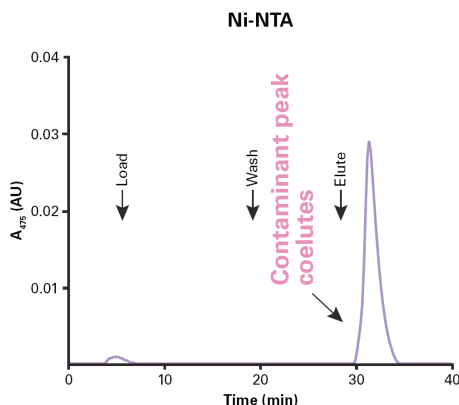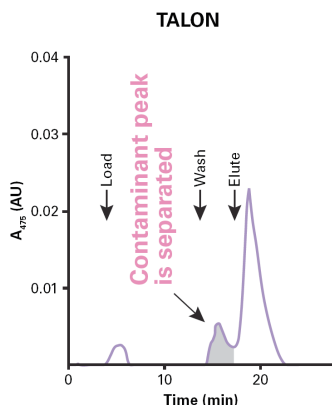IMAC purification—TALON vs. Ni-NTA. 6xHN-AcGFP1 was purified from Sf21 cells using TALON or Ni-NTA. The chromatogram for each column is shown. The absorbance at 475 nm indicates the amount of target protein (AcGFP1) present in each fraction.

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
- Protein labeling
- Immunization to produce antibodies
- Animal Studies


Table I: Reagents Compatible with His60 Ni Superflow Resin

Denaturants	Detergents	Reducing Agents	Others	Salts
6 M Gu-HCl 6 M Urea	2% Triton X-100 2% Tween 20 1% CHAPS	20 mM βME 10 mM DTT	20% glycerol 20% ethanol	4 M MgCl ₂ 5 mM CaCl ₂

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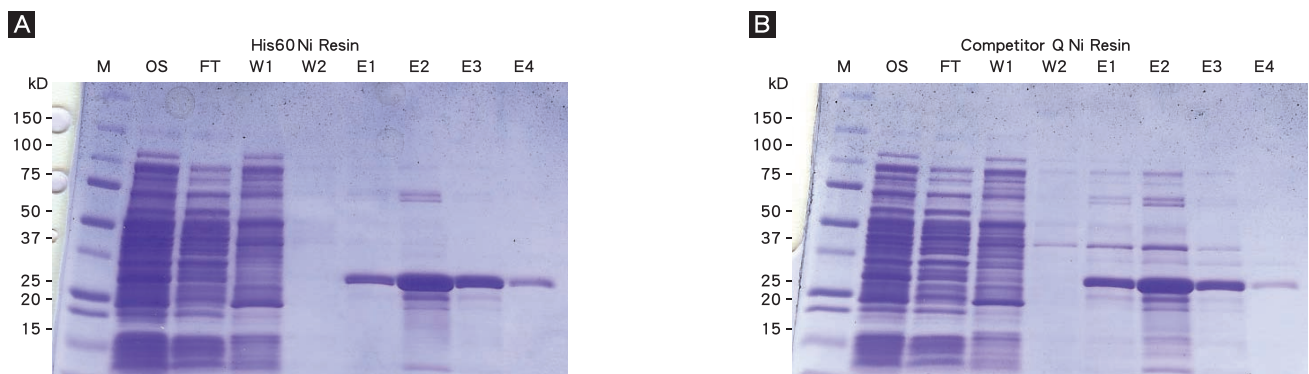
Superior yields and better purity were obtained when using His60 Ni Superflow for purification as compared to Competitor Q's Ni resin. Clontech's pEcoli Linear Expression System was used to express 6xHN-AcGFP1 in *E. coli*. Equivalent amounts of the same sample were used for comparing His60 Ni Superflow Resin purification (Panel A) with Competitor Q Ni Resin purification (Panel B). Lane M: Molecular weight marker. Lane OS: Original sample. Lane FT: Flowthrough. Lanes W1 and W2: Wash. Lane E1-E4: Eluted 6xHN-AcGFP1 fractions.

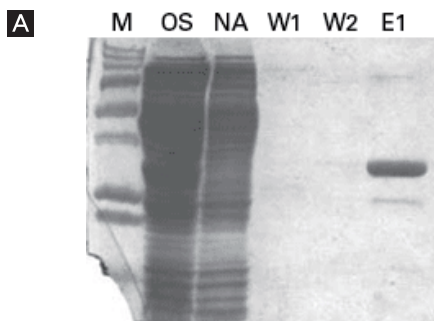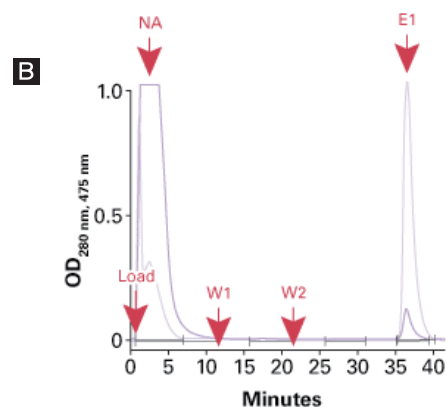A. SDS-PAGE gel analysis of eluate fractions
M: Marker proteins OS: Original sample NA: Nonadsorbed fraction W1: Wash 1 (wash buffer containing 40 mM imidazole) W2: Wash 2 (wash buffer containing 60 mM imidazole) E1: Elution (elution buffer containing 300 mM imidazole)

B. Absorbance at 280 nm/475 nm: imidazole contributes to A280
Load: Loading buffer containing 20 mM imidazole, pH 7.4 W1: Wash buffer containing 40 mM imidazole W2: Wash buffer containing 60 mM imidazole E1: Elution buffer containing 300 mM imidazole

Step gradient elution* results using His60 Ni Superflow Resin.

* Step gradients are recommended over linear gradients for better resolution and sharper elution peaks when eluting from His60 Ni Superflow resin.

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