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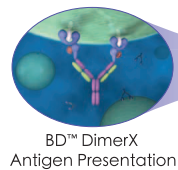
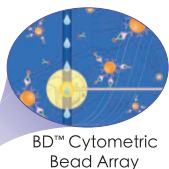
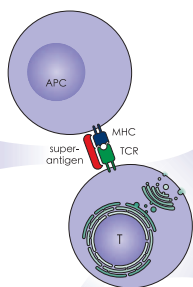
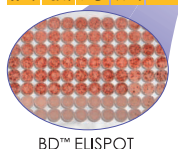
BD™ DimerX Antigen Presentation

BD™ Cytometric Bead Array

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BD™ ELISPOT

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BD™ In Vivo Capture Cytokine Assays

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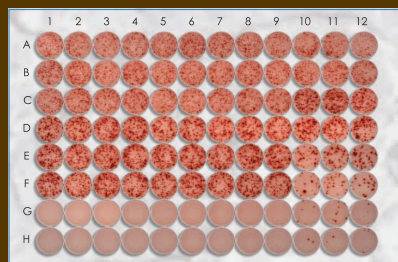
Detection of IFN- $\gamma$ -secreting cells by BD ELISPOT Human IFN- $\gamma$  Kit (Cat. No. 552138). Human PBMCs were stimulated with PMA and ionomycin for 18 hours. (blank controls: G 1-9: cells without stimuli; H 1-9: stimuli without cells)

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**1 Capture Antibody**

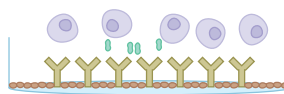
For Sets and Pairs: Coat microwells with anti-cytokine capture antibody.  
For Kits: Go to Step 3; Steps 1 and 2 not necessary.

**2 Blocking**


Block unoccupied sites with protein

**3 Add Cells**


Incubate cells in well with Ag/stimulus etc.

**4 Wash**

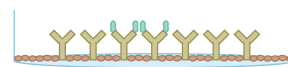
Cells are washed off

**5 Detection Antibody**

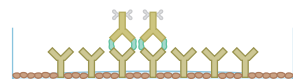
Add Biotinylated anti-cytokine detection antibody

**6 Enzyme-Avidin**

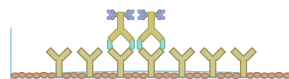Add Avidin-HRP

**7 Develop With Substrate**

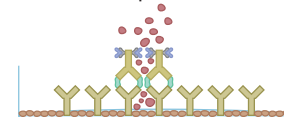Add substrate and monitor formation of colored spots

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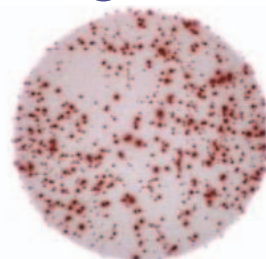
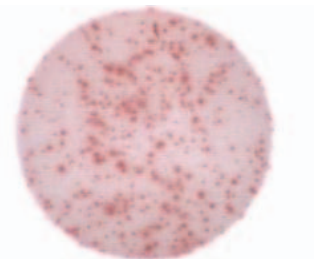
### ELISPOT 包裝

Component	Antibody Pair	Set	Kit
BD ELISPOT plates	未附加	10 uncoated	2 Antibody-coated and blocked
Capture antibody	For 5 plates	For 10 plates	Pre-coated on plates
Detector antibody (biotin-conjugated)	For 5 plates	For 10 plates	For 2 plates
Streptavidin-HRP	需另外購買，貨號：557630	For 10 plates	For 2 plates
AEC chromogen	需另外購買，貨號：551951	需另外購買，貨號：551951	For 2 plates
Wash concentrate, PBS concentrate, assay diluent	未附加	未附加	For 2 plates

\* ELISPOT plates 建議搭配 Millipore MSIPS4510 ( MultiScreenHTS-IP Filter Plate ) for ELISPOT Assay

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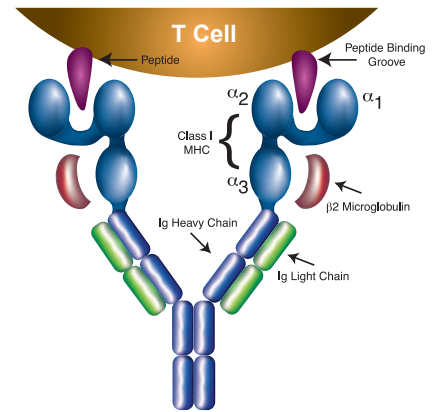
Description	Size	Cat. No.	Description	Size	Cat. No.
Human IL-2 ELISPOT Pair	5 plates	551884	Mouse IL-2 ELISPOT Pair	5 plates	551876
Human IL-2 ELISPOT Set	10 plates	551282	Mouse IL-2 ELISPOT Set	10 plates	551076
Human IL-4 ELISPOT Pair	5 plates	551885	Mouse IL-4 ELISPOT Pair	5 plates	551878
Human IL-4 ELISPOT Set	10 plates	551084	Mouse IL-4 ELISPOT Set	10 plates	551017
Human IL-5 ELISPOT Pair	5 plates	551886	Mouse IL-5 ELISPOT Pair	5 plates	551880
Human IL-5 ELISPOT Set	10 plates	551085	Mouse IL-5 ELISPOT Set	10 plates	551075
Human IL-10 ELISPOT Pair	5 plates	551883	Mouse IL-6 ELISPOT Set	10 plates	552567
Human IL-10 ELISPOT Set	10 plates	551018	Mouse IL-10 ELISPOT Pair	5 plates	557982
Human IL-12p70 ELISPOT Set	10 plates	552574	Mouse IFN- $\gamma$ ELISPOT Pair	5 plates	551881
Human IFN- $\gamma$ ELISPOT Pair	5 plates	551873	Mouse IFN- $\gamma$ ELISPOT Set	10 plates	551083
Human IFN- $\gamma$ ELISPOT Set	10 plates	551849	Mouse IFN- $\gamma$ ELISPOT Kit	2 plates	552569
Human IFN- $\gamma$ ELISPOT Kit	2 plates	552138	Mouse TNF ELISPOT Pair	5 plates	551875
Human TNF ELISPOT Pair	5 plates	551882	Mouse TNF ELISPOT Set	10 plates	551491
Human TNF ELISPOT Set	10 plates	551446	Rat IFN- $\gamma$ ELISPOT Set	10 plates	557819
Human Granzyme B ELISPOT Set	10 plates	552572			

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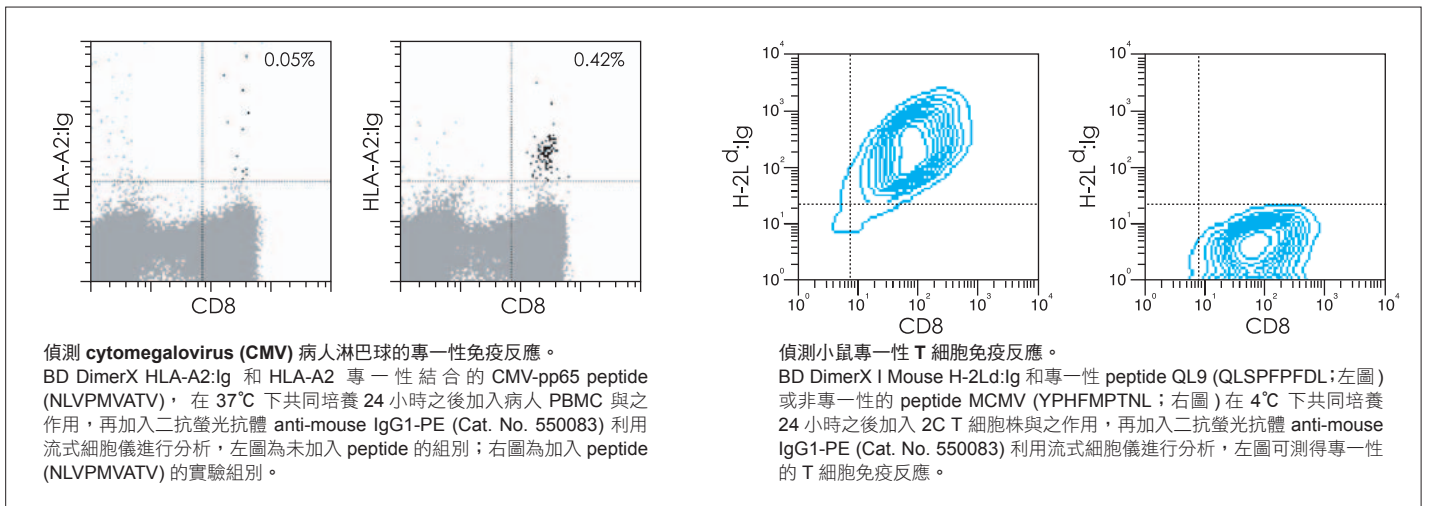
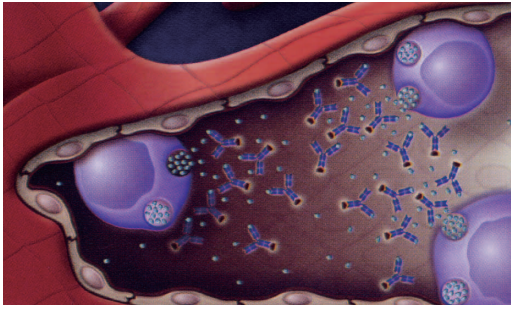
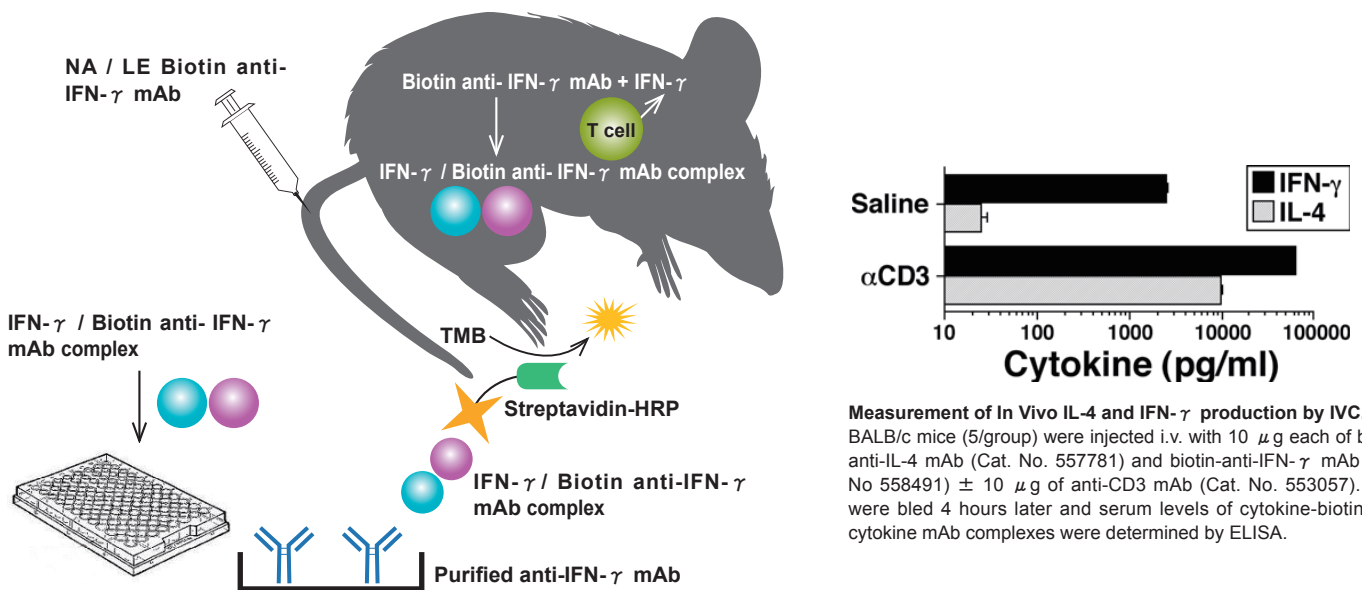
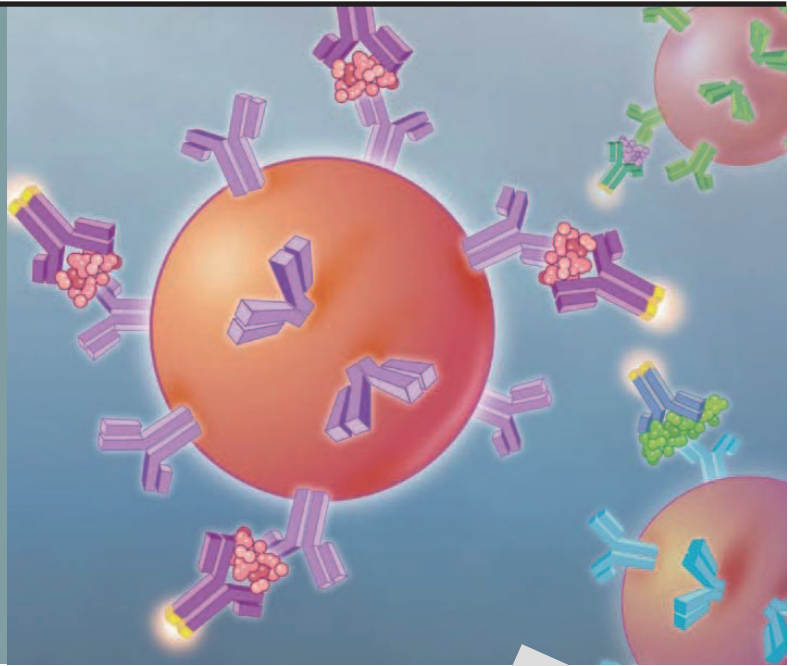
Assay Components :

- *In Vivo* Capture monoclonal antibody (biotinylated, BD NA/LE™)
- Recombinant cytokine standard
- Streptavidin-HRP
- Purified capture monoclonal antibody for plate coating

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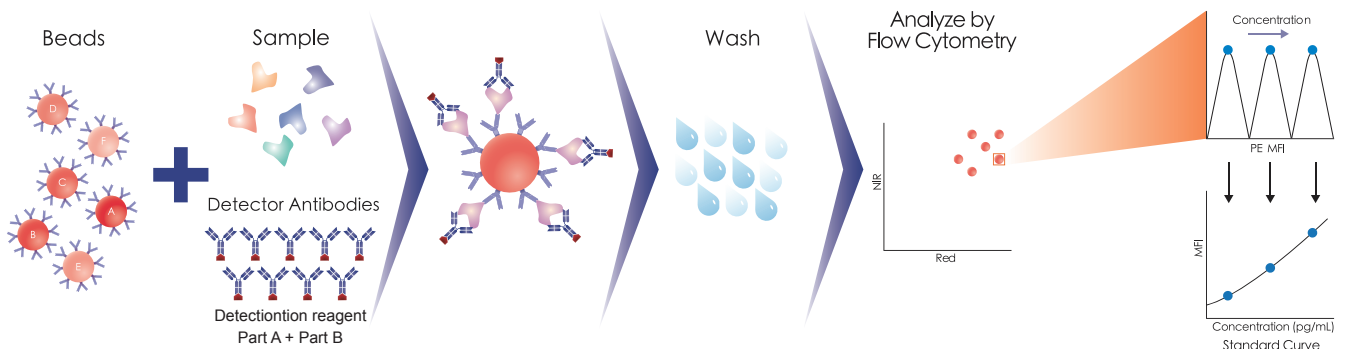
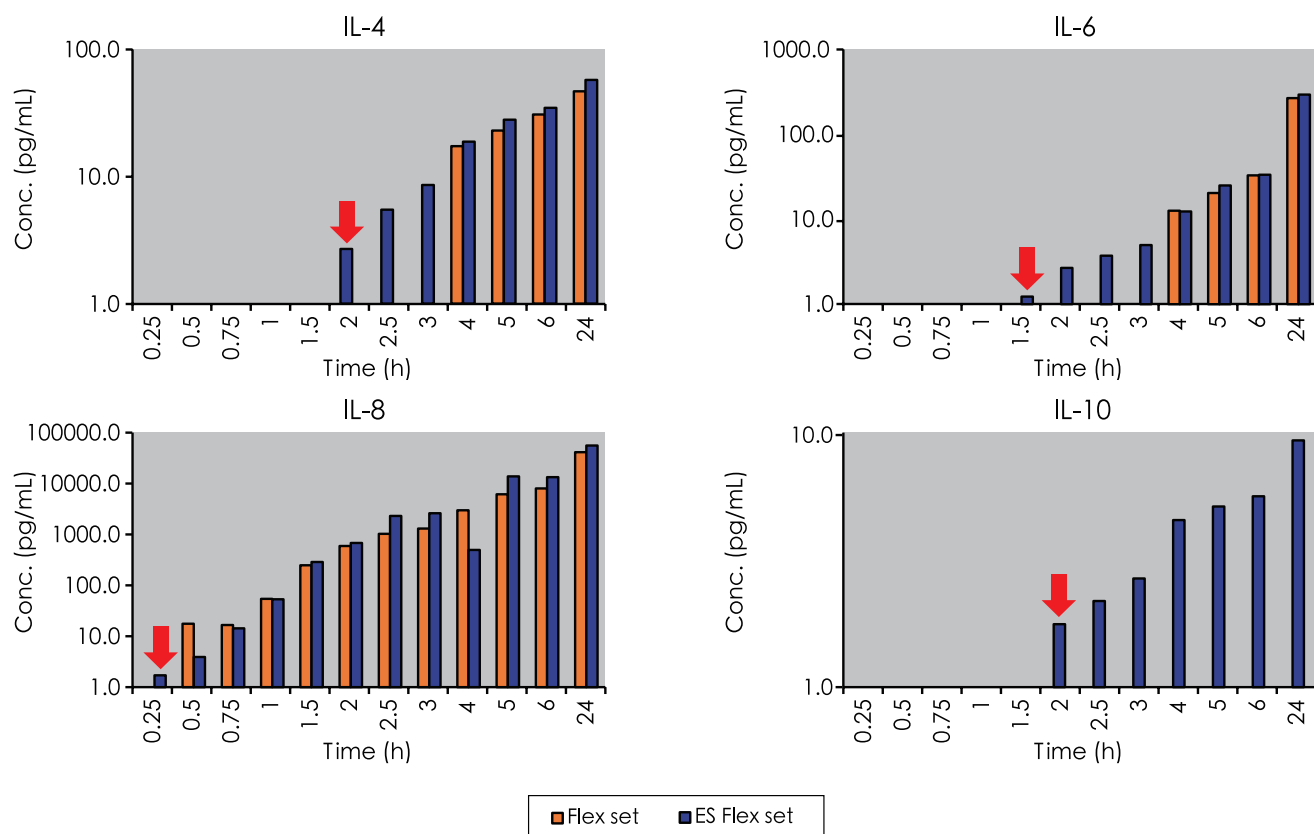
## Time course of human PBMCs stimulated with PMA and Calcium Ionophore.

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