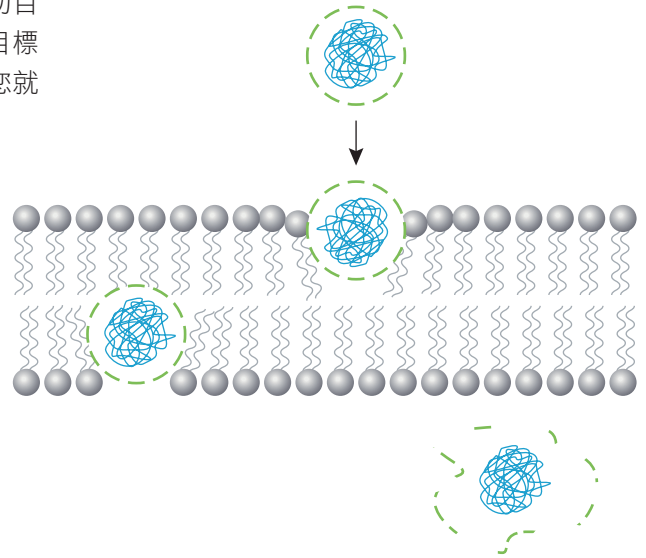
# Xfect™ Protein Transfection Reagent

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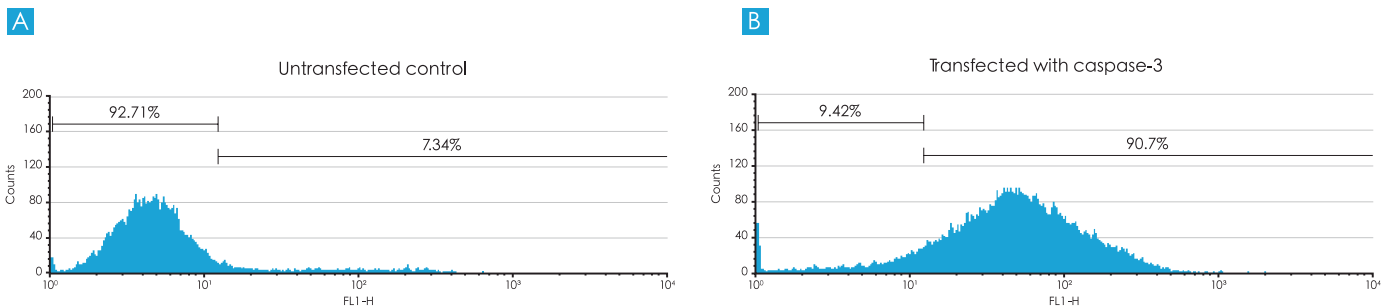
**Xfect Protein Transfection Reagent**  
(cell-penetrating peptide)

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
Complete induction of an early apoptosis event within 2 hours via protein transfection of active caspase-3. Apoptosis is commonly detected using annexin V-FITC staining (see Clontech's ApoAlert™ Annexin V-FITC Apoptosis Kit, Cat. No. 630109), which detects translocation of phosphatidylserine from the inner (cytoplasmic) leaflet of the plasma membrane to the outer (cell surface) leaflet soon after the induction of apoptosis. Annexin V-FITC staining is detected by flow cytometric measurement of increased fluorescence intensity. We transfected human recombinant caspase-3 using Xfect Protein Transfection Reagent and detected completion of apoptosis after just 2 hr, demonstrating that protein transfection using Xfect is both fast and delivers active protein (Panel B). In comparison, other methods for inducing apoptosis require 12 hr for completion of Annexin V staining (1).

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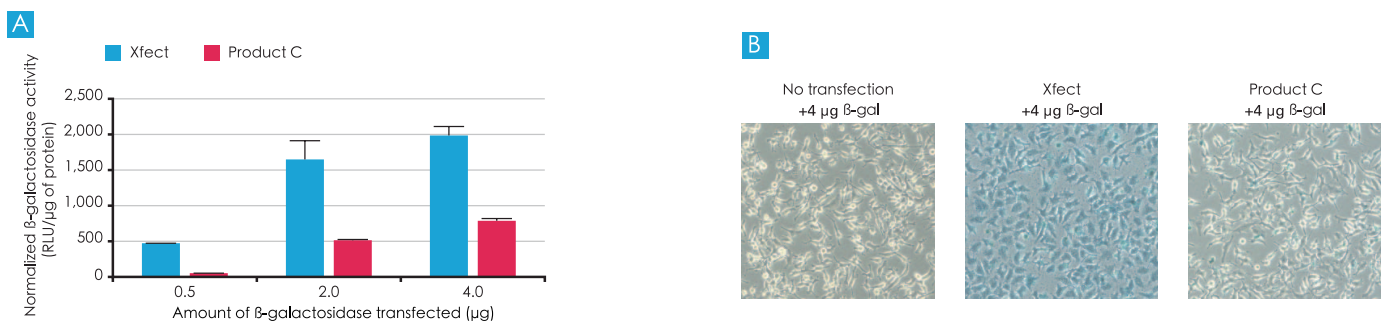Xfect Protein Transfection Reagent delivers up to 8X more active β-galactosidase into HeLa cells than the leading competitor, Product C. Panel A. Xfect Protein Transfection Reagent or Product C was used to transfect Jurkat cells with various amount of β-galactosidase (b-gal) according to the manufacturers' recommended protocol. One hour post-transfection, the cells were assayed for b-gal activity using Clontech's Luminescent β-galactosidase Detection Kit II (Cat. No. 631712). Xfect Protein Transfection Reagent displayed a far higher signal for b-gal than did Product C. Panel B. HeLa cells transfected with 4 μg of b-gal demonstrated higher efficiency and a higher amount of b-gal protein per cell when transfection was performed with Xfect Protein Transfection Reagent (middle panel) than with Product C (righthand panel).