

# The CellASIC® ONIX2 Microfluidic System

Precision control of your cell culture environment  
for advanced live cell imaging & microscopy



The life science business of Merck  
operates as MilliporeSigma in the  
U.S. and Canada.

**MERCK**

# Leading the way in science and technology.

Merck KGaA, Darmstadt, Germany is a leading science and technology company in healthcare, life science and performance materials. Around 50,000 employees work to further develop technologies that improve and enhance life—from biopharmaceutical therapies to treat cancer or multiple sclerosis, to cutting-edge systems for scientific research and production, to liquid crystals for smartphones and LCD televisions. Founded in 1668, we are the world’s oldest pharmaceutical and chemical company.

Our Life Science team is dedicated to providing scientists and engineers with best-in-class technologies, lab materials and services with the intention of making research and production simpler, faster and more successful.

Our solutions enable scientists to spend more time advancing the promise of science through technologies that help detect the previously undetectable, and products that make it possible to monitor live cells in intricate detail.



Cleanroom for CellASIC® ONIX2 Microfluidic System, Hayward, California, USA.

# Quality, reliability and a growing array of publications.

## Introducing the CellASIC® ONIX2 Microfluidic System

With a history of success, dedication to innovation and expertise in life science, Merck KGaA, Darmstadt, Germany, is committed to providing the highest quality, most reliable and relevant products for life science research and production.

The CellASIC® ONIX2 System incorporates extraordinary improvements to traditional live cell observation and experimentation, using advanced microfluidics technology with computer automation to manipulate the culture environment under controlled conditions while allowing you to monitor the cells as they react in real time.

This second generation CellASIC® ONIX2 Microfluidic System is backed by a strong history of distinctive publications, stringent, globally certified manufacturing and quality processes, and a multinational organization committed to delivering the utmost in local service and support.



# Life Is Dynamic.

## Why study living processes by killing your cells?


Scientists recognize that live cell analysis yields necessary and unique insight into living processes that are rapidly changing. While collecting data from a moment in time with classical endpoint analyses has value, the increasingly common practice of observing and measuring cellular changes over extended periods of time reveals more about the true behavior of biological systems. The challenge has been in creating a dynamic cell culture environment that is controllable, manipulatable and reproducible.

“Microfluidic technology is changing the way we perform live cell analysis. The CellASIC® ONIX Microfluidic System is **amazingly easy to use yet highly adaptable**, which greatly reduces the barrier to adopting this emerging technology.”


— Dr. Jintao Liu, Ph.D., San Diego, CA.

## Live cell imaging to transform your research


Move beyond the status quo and gain unprecedented insight into live cells using advanced microfluidics technology. With the CellASIC® ONIX2 Microfluidic System, you too can easily and efficiently create a stable, reproducible environment for imaging. See how it can transform your research.

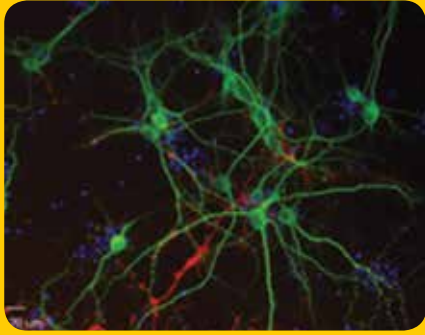

Culturing primary rat cortical neurons (GFP) cultured in the CellASIC® ONIX M04S mammalian microfluidic plate.

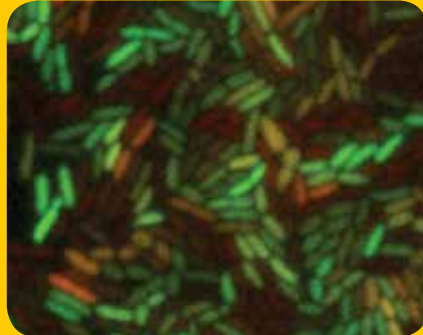

Measuring a gene circuit in *E. coli* in a time-lapse experiment with the CellASIC® ONIX B04 bacterial microfluidic plate. Images were acquired at 100x magnification.

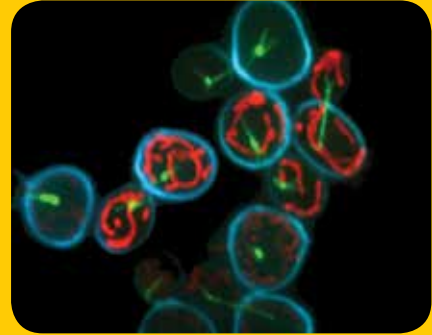

*S. cerevisiae* yeast cells with labeled mitochondria (red), cell wall (cyan), and tubulin (green), imaged in CellASIC® ONIX yeast haploid plate.

Image credit: Maja Bialecka-Fornal, University of California, Irvine, Rafelski Lab.

# The CellASIC® ONIX2 Microfluidic System

Precision control of your cell culture environment for advanced live cell imaging and microscopy.

The ability to grow, observe and manipulate complex cultures requires precision control over the cell culture environment. The second generation CellASIC® ONIX2 Microfluidic System is a refined, yet powerful automated platform for precise manipulation of multiple key cell culture parameters, enabling measurement of cellular responses to pre-programmed media, temperature and gas environment changes.

The CellASIC® ONIX2 Microfluidic System uses high-quality, optically clear microfluidic plates and intuitive software, while integrating with a broad range of inverted microscopes to allow continuous, high magnification observation of live cells — as they react to their environment in time.



1

### Controller System

Small footprint, integrated microincubator controller maintains fluid movement, reagent additions, temperature and gas conditions.

2

### Microfluidic Plates

Application-specific plates bring new cell culture capabilities for live cell imaging.

3

### Manifold


Low-profile, high-quality, high optical clarity culture chambers and manifold mount easily on an inverted microscope—without the need for a bulky environmental chamber.

4

### Software

Intuitive software enables quick and easy setup of detailed protocols for truly automated hands-free cell culture.



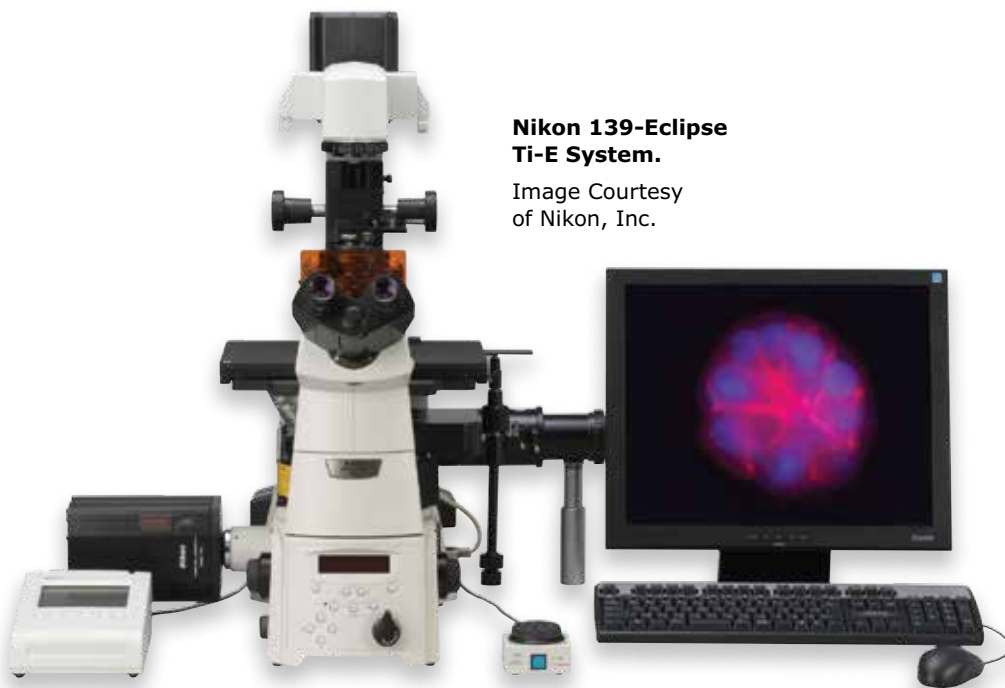

**Nikon 139-Eclipse  
Ti-E System.**

Image Courtesy  
of Nikon, Inc.

## The boost your microscope needs for advanced live cell imaging.

The CellASIC® ONIX2 System turns your microscope into a powerful live cell culture and imaging system that works in conjunction with your capture and analysis software.

You've invested in a powerful microscope; don't limit its use to observing live cells in a static fashion. Enhance your return on investment and obtain more meaningful data using the CellASIC® ONIX2 Microfluidic System as your live cell imaging platform. The low-profile manifold and plate assembly are easily positioned and removed from the microscope stage, providing maximum ease of use and flexibility.

"The key advantage of live cell imaging is that you get to actually watch things unfold before your eyes. It's very powerful to see, visually, what the cells are doing and how they change in space and as a function of time, because biology is a dynamic process...  
**and watching the cells do what they do is something that really appeals to us and is driving our science."**

—Dr. Gurol Suel, San Diego, CA



# There is simply no better way to conduct live cell analysis.

Watch cells change through time and space—all with the precision of the CellASIC® ONIX2 System. Automatically control flow rates, gas and temperature shifts, standing gradients, nutrient/drug additions and media changes. With uninterrupted high-resolution microscopic culture observation and truly consistent, controlled cell culture, you'll answer the questions that set your research apart.

## Eight critical cell culture parameters can be controlled by the CellASIC® ONIX2 Microfluidic System:

Parameter	If too low, can cause:	If too high, can cause:
Temperature	Decreased cell response	Increased respiration/protein damage
Oxygen Level	Decreased pH/increased glycolysis	Increased ROS, membrane damage
Growth Factors	Increased apoptosis/decreased protein synthesis	Increased angiogenesis and cell division
Humidity	Increased osmolarity/cell metabolism/oxidative stress	Could damage imaging equipment
pH	Protein and membrane denaturation	Increased alkalosis and dehydration
Osmolality	Decreased cell division/increased autophagic proteolysis and cell rupture	Increased oxidative stress, DNA breakage and nutrient digestion
Glucose	Decreased autophagy and metabolism	Increased apoptosis and ROS
ECM and Adhesion	Decreased angiogenesis/aberrant differentiation	Increased cell adhesion, chemotaxis, proliferation



Small footprint, integrated microincubator controller maintains fluid movement, reagent additions, temperature and gas mixture.


## Application-specific microfluidic plates for optimized, bio-inspired cell culture.

The key to microenvironment control using the CellASIC® ONIX2 System is in coupling precision-manufactured microfluidic cell culture plates with an automated perfusion system to provide a controlled environment optimized for high-resolution live cell imaging.

### Optimized, bio-inspired cell culture

Different cells need different environments. CellASIC® ONIX microfluidic plates are designed to optimize cell health during dynamic live cell experiments.

Various application specific plate designs give you the flexibility to probe the questions that interest you the most.

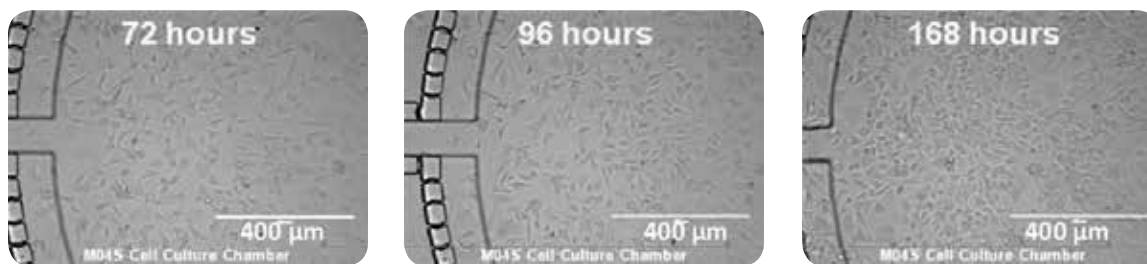


**Healthy long-term cultures outside the incubator.**

NIH 3T3 cells were cultured in the CellASIC® ONIX Microfluidic System (M04S plate) with continuous perfusion and monitored using bright field microscopy for 168 hours.


### Advantages of the CellASIC® ONIX2 microfluidic plates:

- Rapid perturbations of environmental parameters due to the tiny microfluidic volumes involved
- Perfusion microbarriers allow continuous reagent transport without shearing stress
- High-resolution imaging through optical glass bottom and chamber structures that keep cells in a single focal plane
- Ability to run multiple independent experiments simultaneously
- Continuous access of nutrients and ongoing removal of waste, promoting optimal cell health

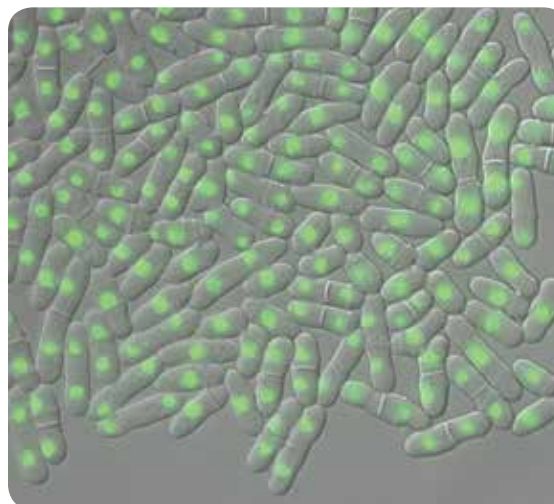

**Fission yeast, cultured and imaged in the CellASIC® Y04D microfluidic plate.**

Image courtesy of Prof. Hironori Niki, National Institute of Genetics, Japan.

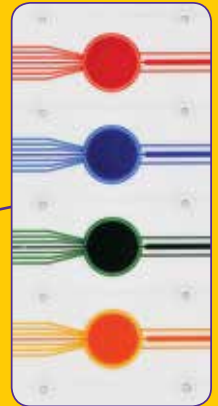

Application-specific plates bring new cell culture capabilities for live cell imaging.


**Merck's strength in engineering and quality systems are ideally suited to meet the challenges of consistent and scalable microfluidic plate manufacturing:**

- Patented microfluidic manufacturing method
- Multi-layer molding capability from 0.5 – 150 microns
- Commitment to continuous improvement of product quality

## Low-profile manifold replaces bulky environmental chambers.

A challenge to high resolution, long-term cell imaging lies in the need to control the culture environment without impeding optical access to the cells. This is complicated by the proximity of microscope elements (objective lens, condenser, stage), and the need for an unobstructed light path through the sample that does not hinder stage or objective movement. The miniaturization provided by the CellASIC® ONIX2 Microfluidic System is not only valuable for easily controlling temperature and fluids, but also maintains a small footprint for use directly on the microscope stage for optimal clarity and ease of use.

A low-profile manifold connects the control system with the microfluidic plates. Various manifold layouts are available to address different plate formats and application types.

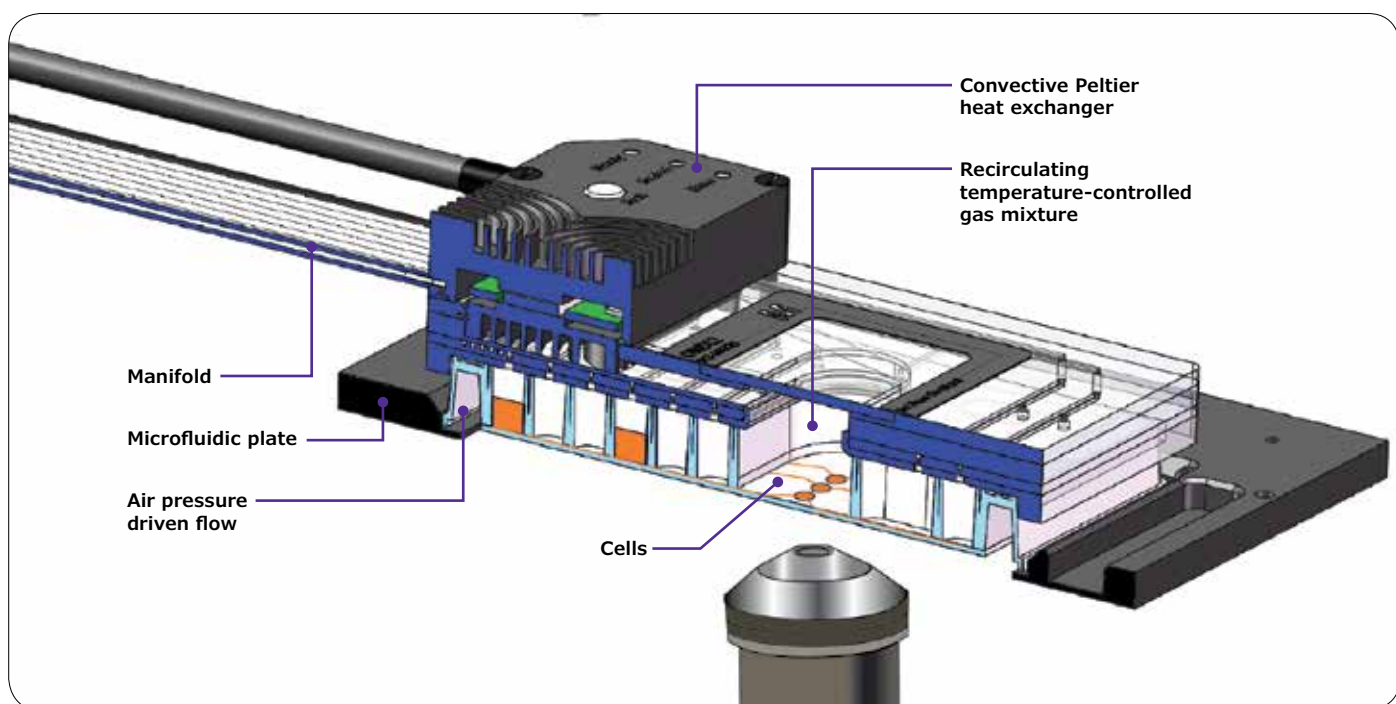

Low-profile, high-quality, high optical clarity culture chambers and manifold mount easily on an inverted microscope—without the need for a bulky environmental chamber.

“What the CellASIC® system lets us do is very rapidly turn on and off conditions or insults while following single cells.”

— Dr. Ethan Garner, Cambridge, MA

#### **Manifold advantages:**

- Manifold seals to the microfluidic plate via “one-touch” vacuum mechanism
- Pressure-driven flow of liquids to the cell chamber provides high precision, even at very small volumes
- No contact between the flow system and the biological solutions (on the plate) prevents contamination
- Low-profile manifold is easy to adapt to any inverted microscope stage

## Intuitive software adds the power of automation to your experiment.

Set your parameters, time your changes and the system does the rest. Simple and straightforward, the intuitive interface makes it possible for experts or novice users to get started quickly and easily, while ensuring consistency throughout the experiment.

### Advanced software provides:

- Clearly defined protocols with all experimental parameters saved in one file
- Convenience and reproducibility. To repeat an experiment, just put in a new plate, and click “run”
- Consistency and reliability. Better portability and control of software protocols means more consistent experimentation
- Error checking, self test and data logging all make for better troubleshooting and serviceability
- The ability to use with your microscopy software. Synchronize imaging with cell experiment control and connect your experiment data with your image data



Intuitive software enables quick and easy setup of detailed protocols for truly automated hands-free cell culture.




“We’ve been able to quickly and easily perform novel and technologically demanding experiments without any prior microfluidic experience. I’ve been able to focus on the fundamental biological questions while letting CellASIC® provide me with the tools I need to answer them.”

— Maheshri Lab, Cambridge, MA

# Discover new insights into dynamic cellular processes.

The CellASIC® ONIX2 Microfluidic System provides detailed information about live cell processes with ultra-high-quality images and complete environment control.

## Dynamic assays on adherent and non-adherent cells

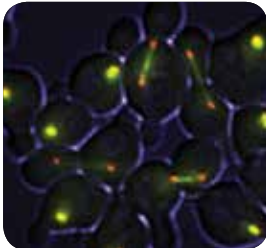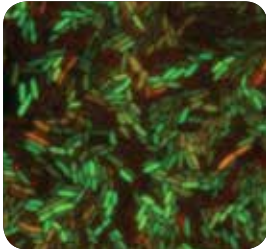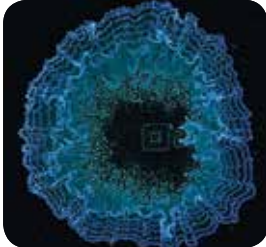CellASIC® ONIX2 Advantage	Application		
High-resolution imaging in single focal plane with controlled nutrient flow for tracking single cells over time	Yeast single cell response		<i>S. cerevisiae</i> cells expressing GFP-tubulin and SPC42- mCherry during alpha-factor exposure and arrest. Images were acquired at 60x magnification. Courtesy of S. Lacefield, U. Indiana.
Measure multi-generational responses to live bacteria while maintaining cells in a single focal plane for days	Bacterial single cell response		A gene circuit in <i>E. coli</i> was induced and visualized for a time-lapse experiment in the CellASIC® ONIX B04 microfluidic plate. Images were acquired at 100x magnification.
Continuous monitoring of community dynamics and precise control of growth environment	Bacterial biofilm dynamics		Time-lapse, composite image of bacterial biofilm growth. Courtesy of the Suel lab, UCSD.






**The CellASIC® ONIX2 Microfluidic System works with your inverted microscope to enable long-term cell perfusion experiments and dynamic time-lapse analysis.**

(Shown with Lionheart™ FX automated Live Cell Imager from BioTek® Instruments)






# Control cell culture conditions to induce and observe reactions in real time.

## Rapid gas and media switching capabilities

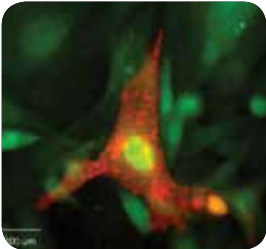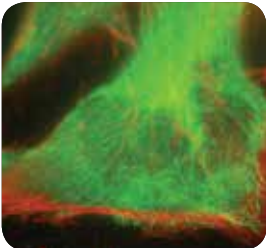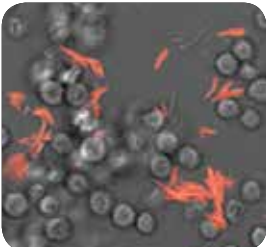CellASIC® ONIX2 Advantage	Application		
Fast and precise control over culture conditions to induce hypoxic, starvation or toxic microenvironments	Cell response to hypoxic environments		LAMP1-RFP/ LC3-GFP CHO reporter cells cultured on the CellASIC® ONIX system showing autophagosomes (green) and lysosomes (red) during a hypoxia-induced autophagy assay over 24 hours.
Microfluidic control of nutrient additions during continuous observation	Cell response to changing media conditions, drugs and other stimulants		Long-term live cell microscopy of cellular cytoskeletal changes in HeLa cells with precise microenvironment control. Cells were stained for tubulin (green) and actin (red) using "in-plate" immunostaining with multi-solution, automated washing and exposure programs, in the CellASIC® ONIX M04S Microfluidic Plate. Image was acquired at 100x magnification.
High-resolution imaging of dynamic cellular interactions	Host-pathogen interactions		Host-pathogen assay monitoring of M. tuberculosis-RFP infection in macrophages.

# Enhance live cell imaging with stable, optimized culture conditions.

## Maintenance of constant environment for optimal cell culture

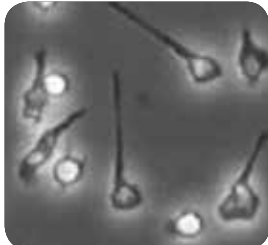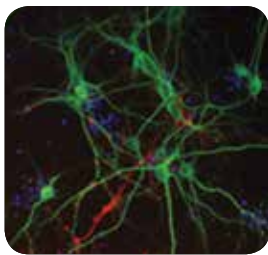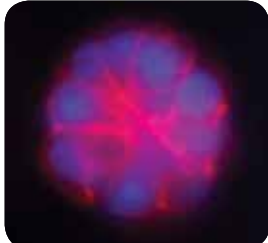CellASIC® ONIX2 Advantage	Application	
<p>Microfluidic control creates stable standing chemical gradients</p>	<p>Chemotaxis/migration in response to chemogradient</p>	 <p>Chemotaxis/migration in response to chemogradient HL-60 neutrophil migration in response to a chemokine. This frame from a live cell analysis video shows cells concentrating toward the chemokine in one chamber of a CellASIC® ONIX M04G Microfluidic Gradient Plate. Courtesy of Jason Park, Wendell Lim Lab, UCSF.</p>
<p>Microfluidics and computer control allow quick adjustments to microenvironment parameters and media changes, all while recording cell behavior</p>	<p>Setting optimal culture parameters for difficult cell types or long-term experiments</p>	 <p>Primary rat cortical neurons cultured in the CellASIC® ONIX M04S Microfluidic Plate to Day 15 and immunostained in-plate for MAP2 (Green GFP, neurons) and GFAP (Red RFP, astrocytes; 40x).</p>
<p>Control and perfusion dynamics create more <i>in vivo</i>-like environment for complex cultures</p>	<p>Microscopy of 3D cell culture</p>	 <p>Observation of multi-day morphology changes of 3D cancer spheroids cultured in extracellular matrix. MCF-10A breast cancer cells were suspended in Matrigel® substrate and grown in the CellASIC® ONIX M04S Microfluidic Plate. Cells were stained for actin (red) and nuclei (blue). Image was acquired at 40x magnification.</p>



“Since I aim to quantify mitochondrial morphology, I require constant, stable imaging conditions that maintain the health of the cells, which the CellASIC® ONIX System does very well.”

- Marshall Lab, San Francisco, CA

## A Growing Array of Distinguished Publications

The CellASIC® ONIX2 Microfluidic System is the next generation of a well-established and highly published technology. Here is a sampling of recent publications.

Find more by visiting our website at: [merckmillipore.com/cellasic-publications](http://merckmillipore.com/cellasic-publications)

### Yeast

Kabeche R, Howard L, Moseley JB. Eisosomes provide membrane reservoirs for rapid expansion of the yeast plasma membrane. *J. Cell Sci.*, Nov 2015; 128: 4057 - 4062.

<http://www.ncbi.nlm.nih.gov/pubmed/26403204>

Peroza EA, Ewald JC, Parakkal G, Skotheim JM, Zamboni N; A genetically encoded FRET sensor for monitoring in vivo trehalose-6-phosphate dynamics; *Analytical Biochemistry* 2015, Apr 1; 474:1-7. doi:10.1016/j.ab.2014.12.019

<http://www.ncbi.nlm.nih.gov/pubmed/25582303>

Kabeche R, Madrid M, Cansado J, Moseley JB. Eisosomes Regulate Phosphatidylinositol 4,5-Bisphosphate (PI(4,5)P2) Cortical Clusters and Mitogen-activated Protein (MAP) Kinase Signaling upon Osmotic Stress. *J. Biol. Chem.*, Oct 2015; 290: 25960 - 25973.

<http://www.ncbi.nlm.nih.gov/pubmed/26359496>

Mazo-Vargas A, Park H, Aydin M, Buchler NE; Measuring fast gene dynamics in single cells with time-lapse luminescence microscopy; *Mol. Biol. Cell* November 5, 2014 vol. 25 no. 22 3699-3708; doi: 10.1091/mbc.E14-07-1187

<http://www.ncbi.nlm.nih.gov/pubmed/25232010>

Burke TA, Christensen JR, Barone E, Suarez C, Sirotkin V, Kovar DR. Homeostatic actin cytoskeleton networks are regulated by assembly factor competition for monomers. *Curr Biol.* 2014 Mar 3;24(5):579-85.

<http://www.ncbi.nlm.nih.gov/pubmed/24560576>

Meyer RE, Kim S, Obeso D, Straight PD, Winey M, Dawson DS. Mps1 and Ipl1/Aurora B act sequentially to correctly orient chromosomes on the meiotic spindle of budding yeast. *Science.* 2013 Mar 1;339(6123):1071-4.

<http://www.ncbi.nlm.nih.gov/pubmed/23371552>

Rafelski SM, Viana MP, Zhang Y, Chan YH, Thorn KS, Yam P, Fung JC, Li H, Costa L da F, Marshall WF. Mitochondrial network size scaling in budding yeast. *Science.* 2012 Nov 9;338(6108):822-4.

<http://www.ncbi.nlm.nih.gov/pubmed/23139336>

Kraft C, Kijanska M, Kalie E, Siergiejuk E, Lee SS, Semplicio G, Stoffel I, Brezovich A, Verma M, Hansmann I, Ammerer G, Hofmann K, Tooze S, Peter M. Binding of the Atg1/ULK1 kinase to the ubiquitin-like protein Atg8 regulates autophagy. *EMBO J.* 2012 Sep 12;31(18):3691-703.

<http://www.ncbi.nlm.nih.gov/pubmed/22885598>

Kono K, Saeki Y, Yoshida S, Tanaka K, Pellman D. Proteasomal degradation resolves competition between cell polarization and cellular wound healing. *Cell.* 2012 Jul 6;150(1):151-64.

<http://www.ncbi.nlm.nih.gov/pubmed/22727045>

Sanchez-Diaz A, Nkosi PJ, Murray S, Labib K. The Mitotic Exit Network and Cdc14 phosphatase initiate cytokinesis by counteracting CDK phosphorylations and blocking polarised growth. *EMBO J.* 2012 Aug 29;31(17):3620-34.

<http://www.ncbi.nlm.nih.gov/pubmed/22872148>

Wei P, Wong WW, Park JS, Corcoran EE, Peisajovich SG, Onuffer JJ, Weiss A, Lim WA. Bacterial virulence proteins as tools to rewire kinase pathways in yeast and immune cells. *Nature.* 2012 Aug 16;488(7411):384-8.

<http://www.ncbi.nlm.nih.gov/pubmed/22820255>

Bermejo C, Haerizadeh F, Takanaga H, Chermak D, Frommer WB. Optical sensors for measuring dynamic changes of cytosolic metabolite levels in yeast. *Nat Protoc.* 2011 Oct 27;6(11):1806-17.

<http://www.ncbi.nlm.nih.gov/pubmed/22036883>

Eser U, Falleur-Fettig M, Johnson A, Skotheim JM. Commitment to a cellular transition precedes genome-wide transcriptional change. *Mol Cell.* 2011 Aug 19;43(4):515-27.

<http://www.ncbi.nlm.nih.gov/pubmed/21855792>

Dechant R, Binda M, Lee SS, Pelet S, Winderickx J, Peter M. Cytosolic pH is a second messenger for glucose and regulates the PKA pathway through V-ATPase. *EMBO J.* 2010 Aug 4;29(15):2515-26.

<http://www.ncbi.nlm.nih.gov/pubmed/20581803>

Manzoni R, Montani F, Visintin C, Caudron F, Ciliberto A, Visintin R. Oscillations in Cdc14 release and sequestration reveal a circuit underlying mitotic exit. *J Cell Biol.* 2010 Jul 26;190(2):209-22.

<http://www.ncbi.nlm.nih.gov/pubmed/20660629>

Furuya K, Niki H. The DNA damage checkpoint regulates a transition between yeast and hyphal growth in *Schizosaccharomyces japonicus*. *Mol Cell Biol.* 2010 Jun;30(12):2909-17. doi: 10.1128/MCB.00049-10.

<http://www.ncbi.nlm.nih.gov/pubmed/20368354>

Octavio LM, Gedeon K, Maheshri N. Epigenetic and conventional regulation is distributed among activators of FLO11 allowing tuning of population-level heterogeneity in its expression. *PLoS Genet.* 2009 Oct;5(10):e1000673.

<http://www.ncbi.nlm.nih.gov/pubmed/19209350>

## Bacteria

Sutterlin HA, Shi H, May KL, Miguel A, Khare S, Huang KC, and Silhavy TJ. Disruption of lipid homeostasis in the Gram-negative cell envelope activates a novel cell death pathway. PNAS. 2016 Feb; 10.1073/pnas.1601375113.

<http://www.ncbi.nlm.nih.gov/pubmed/26929379>

Prindle A, Liu J, Asally M, Ly S, Garcia-Ojalvo J, Süel GM. Ion channels enable electrical communication in bacterial communities. Nature. 2015 Nov 5;527(7576):59-63.

<http://www.ncbi.nlm.nih.gov/pubmed/26503040>

Grangeon R, Zupan JR, Anderson-Furgeson J, and Zambryski PC. PopZ identifies the new pole, and PodJ identifies the old pole during polar growth in *Agrobacterium tumefaciens*. PNAS. 2015 Sep; 112:11666 - 11671.

<http://www.ncbi.nlm.nih.gov/pubmed/26324921>

Liu J, Prindle A, Humphries J, Gabalda-Sagarra M, Asally M, Lee DD, Ly S, Garcia-Ojalvo J, Süel GM. Metabolic co-dependence gives rise to collective oscillations within biofilms. Nature 2015 July 523:550-554.

<http://www.ncbi.nlm.nih.gov/pubmed/26200335>

Sieger B, Schubert K, Donovan C, Bramkamp M. The lipid II flippase RodA determines morphology and growth in *Corynebacterium glutamicum*. Mol Microbiol. 2013 Dec;90(5):966-82.

<http://www.ncbi.nlm.nih.gov/pubmed/24118443>

Gordon AJ, Satory D, Halliday JA, Herman C. Heritable change caused by transient transcription errors. PLoS Genet. 2013 Jun;9(6):e1003595.

<http://www.ncbi.nlm.nih.gov/pubmed/23825966>

Donovan C, Schauss A, Kramer R, Bramkamp M. Chromosome segregation impacts on cell growth and division site selection in *Corynebacterium glutamicum*. PLOS One, February 2013; 8(2): e55078.

<http://www.ncbi.nlm.nih.gov/pubmed/23405112>

Enrique Rojas, Julie A. Theriot, and Kerwyn Casey Huang. Response of *Escherichia coli* growth rate to osmotic shock. PNAS, May 2014; 111: 7807 - 7812.

<http://www.ncbi.nlm.nih.gov/pubmed/24821776>

Young JW, Locke JC, Elowitz MB. Rate of environmental change determines stress response specificity. Proc Natl Acad Sci U S A. 2013 Mar 5;110(10):4140-5.

<http://www.ncbi.nlm.nih.gov/pubmed/23407164>

## Algae

Ludington WB, Shi LZ, Zhu Q, Berns MW, Marshall WF. Organelle size equalization by a constitutive process. Curr Biol. 2012 Nov 20;22(22):2173-9.

<http://www.ncbi.nlm.nih.gov/pubmed/23084989>

## Mammalian

Changou CA, Chen Y-R, Xing L, Yen Y, Chuang FYS, Cheng RH, Bold RJ, Ann DK, Kung H-J; Arginine starvation-associated atypical cellular death involves mitochondrial dysfunction, nuclear DNA leakage, and chromatin autophagy. PNAS, Sep 2014; 111: 14147 - 14152.

<http://www.ncbi.nlm.nih.gov/pubmed/25122679>

Zambrano S, De Toma I, Piffer A, Bianchi ME, Agresti A; NF- $\kappa$ B oscillations translate into functionally related patterns of gene expression; eLife 2016; 10.7554/eLife .09100.

<http://www.ncbi.nlm.nih.gov/pubmed/26765569>

Park JS, Rhau B, Hermann A, McNally KA, Zhou C, Gong D, Weiner OD, Conklin BR, Onuffer J, Lim WA; Synthetic control of mammalian-cell motility by engineering chemotaxis to an orthogonal bioinert chemical signal. PNAS, Apr 2014; 111: 5896 - 5901.

<http://www.ncbi.nlm.nih.gov/pubmed/24711398>



View the updated list of publications, review protocols and application data and watch video of live cells responding in real time by visiting: [merckmillipore.com/cellasic](http://merckmillipore.com/cellasic)

## Ordering Information

Description	Qty	Catalog No.
<b>CellASIC® ONIX2 Microfluidic System and Manifolds</b>		
CellASIC® ONIX2 Microfluidic System	1	<b>CAX2-S0000</b>
CellASIC® ONIX2 Manifold XT (Temperature Controlled)	1	<b>CAX2-MXT20</b>
CellASIC® ONIX2 Manifold Basic (No Temperature Control)	1	<b>CAX2-MBC20</b>
<b>Microfluidic Plates</b>		
CellASIC® ONIX Plate for Haploid Yeast Cells	5	<b>Y04C-02-5PK</b>
CellASIC® ONIX Plate for Diploid Yeast Cells	5	<b>Y04D-02-5PK</b>
CellASIC® ONIX Plate for Bacteria Cells	5	<b>B04A-03-5PK</b>
CellASIC® ONIX Switching Plate for Mammalian Cells	5	<b>M04S-03-5PK</b>
CellASIC® ONIX Gradient Plate for Mammalian Cells	5	<b>M04G-02-5PK</b>
CellASIC® ONIX Open-Top Plate for Mammalian Cells	5	<b>M04L-03-5PK</b>
CellASIC® ONIX Plate for <i>Chlamydomonas</i> Cells	5	<b>C04A-01-5PK</b>



For a complete list of products and information, visit: [merckmillipore.com/cellasic](https://www.merckmillipore.com/cellasic)



# Technical Specifications

General	
Dimensions	System: 330 mm wide × 306 mm deep × 108 mm high (13 in. × 12 in. × 4.25 in.)
Weight	System: 6.7 kg (14.75 lb)
Power consumption	100–240 VAC, 50/60 Hz 40 W
Cooling mode	Air cooled (natural convection)
Environmental conditions	For indoor laboratory use
Operating temperature	20°C to 30°C
Storage temperature	5°C to 35°C
Stability of sample chamber temperature	± 0.2°C
Rise time (25°C to 37°C)	< 30 minutes
Cooling time (37°C to 25°C)	< 30 minutes
Flow Control	
Number of outputs	8 (each addressable by either of two pressure controllers)
Pressure range	-50 to 70 kPa (-7.25–10.2 psi)
Pressure accuracy	± 1.5 kPa (0.22 psi)
Pressure stabilization time	within ± 5 kPa (0.73 psi) in < 5 seconds
Gas Environment Control	
Control capabilities	Option to electronically select between two premixed gasses and set flow rate to slow or fast
Gas input requirements	Clean, dry, premixed gas mixtures containing air, CO <sub>2</sub> , N <sub>2</sub> , and oxygen (up to 25%), regulated to between 100 kPa and 700 kPa (15 psi and 100 psi)
Gas consumption	Slow flow: 5 mL/min ± 2 mL/min Fast flow: 50 mL/min ± 20 mL/min
Cell culture region gas environment accuracy	For gas flow at 3 mL/min: < 10% deviation from delivered gas concentration For gas flow at 30 mL/min: < 2% deviation from delivered gas concentration
Temperature Control	
Temperature control range	Room temperature to 40°C
Control method	Bi-directional PID
Accuracy of sample chamber temperature	± 1°C (using the CAX2-ACT20 ONIX2 Temperature Calibration Plate can result in accuracy as high as ± 0.2°C)
Computer Requirements	
Operating system	Windows® 7, or Windows® 10 (32 or 64 bit)
Random access memory (RAM)	4 GB or higher
Monitor	1920x1080 resolution or higher
Hard drive	200 MB or higher
Hardware interface	USB 2.0 or higher
Microscope Requirements	
Type	Inverted microscope
Stage weight capacity	For CAX2-MXT20 CellASIC® ONIX2 Manifold XT and filled plate: 325 g (11.1 oz) For CAX2-MBC20 CellASIC® ONIX2 Manifold Basic and filled plate: 150 g (5.3 oz)
Minimum condenser working distance	For CAX2-MXT20 CellASIC® ONIX2 Manifold XT: 28 mm (0.9 in) For CAX2-MBC20 CellASIC® ONIX2 Manifold Basic: 28 mm (0.9 in)
Maximum condenser diameter	For CAX2-MXT20 CellASIC® ONIX2 Manifold XT: 70 mm (2.8 in) For CAX2-MBC20 CellASIC® ONIX2 Manifold Basic: no limit

# Cultivate Consistency

Create a consistent cell culture environment with our comprehensive portfolio of reagents and cultureware. From cell preparation and growth to investigation and analysis, our products provide the quality and reliability to get the most out of your time and research. Explore our full line of cell culture solutions to further your cellular research with confidence.

Learn more at: [sigmaaldrich.com/cellculture](https://sigmaaldrich.com/cellculture)

**Streamline  
your cell culture  
experiments and get  
the most from your  
CellASIC® ONIX2  
Microfluidic System.**



## PREPARE

### Type of Cells

- **Cell Lines**  
Cell lines can be cultured indefinitely and are therefore versatile, cost-effective tools for modeling more complex biological systems. Count on the highest quality cell lines through our partnership with the European Collection of Authenticated Cell Cultures (ECACC) delivering the most diverse selection from which to choose.
- **Primary Cells**  
Closely mimic what occurs *in vivo* using our extensive collection of primary cells that have been isolated directly from human or animal tissue.

### Filtration

Sterilize your cell culture media to eliminate contaminants and ensure reproducibility with our complete line of sterile filtration products, including the one and only Stericup® Vacuum Filtration System.

### Cell freezing solutions

Ensure proper preservation of your cells by using our application-tested cryoprotectants and ready-to-use freezing media.

## GROW

### Liquid Media & Supplements

Support the growth of your cells with our comprehensive offering of classical and specialty media. We also carry a wide variety of application-tested supplements and growth factors to meet all your research needs.

### Specialty Culture Inserts & Plates

Promote cell growth with access to media from both the apical and basolateral sides to closely mimic what occurs *in vivo* with Millicell® cell culture inserts and plates. Choose from standing, hanging, or organotypic insert formats, available with a range of membranes to deliver the growth environment that suits your application.

Our MultiScreen® plates (24- or 96-well) have a long history as a reliable tool for the advancement of drug discovery and life science research in multiple applications.

### Corning® Cultureware

Support your cell growth with a comprehensive line of high-quality cell culture and filtration solutions from Corning®, including serological pipettes, cell culture multiple well plates, flasks, dishes, permeable supports and filters.


## ANALYZE

### Scepter™ 2.0 Cell Counter

Simplify your cell counting with an automated, hand-held device that provides reliable cell counts in less than 30 seconds.

### CellASIC® ONIX2 Microfluidic System

The CellASIC® ONIX2 Microfluidic System automates your cell culture. Using high-quality, optically clear microfluidic plates and intuitive software, the system integrates with a broad range of inverted microscopes to allow continuous, high magnification observation of live cells — as they react to their environment in real time.

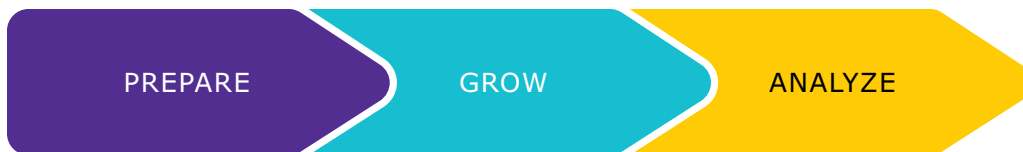

- Cell Lines & Primary Cells
- Filtration
- Cryopreservation

- Liquid Media & Supplements
- Specialty Culture Inserts & Plates
- Corning® Cultureware

- Scepter™ 2.0 Cell Counter
- CellASIC® ONIX2 Microfluidic System

**Merck**  
Streamline your cell culture workflow.

# The CellASIC® ONIX2 Microfluidic System



To request a demonstration, or for more information on the CellASIC® ONIX2 System, please visit: [merckmillipore.com/cellasic](https://www.merckmillipore.com/cellasic)

To learn more about live cell imaging and microfluidics technology, visit our learning center at: [merckmillipore.com/cellasic-livecell](https://www.merckmillipore.com/cellasic-livecell)



**To place an order or receive technical assistance**

In Europe, please call Customer Service:

France: 0825 045 645

Germany: 069 86798021

Italy: 848 845 645

Spain: 901 516 645 Option 1

Switzerland: 0848 645 645

United Kingdom: 0870 900 4645

For other countries across Europe,  
please call: +44 (0) 115 943 0840

Or visit: [www.merckmillipore.com/offices](http://www.merckmillipore.com/offices)

For Technical Service visit:

[www.merckmillipore.com/techservice](http://www.merckmillipore.com/techservice)

[merckgroup.com/life-science](http://merckgroup.com/life-science)

