Liquid Biopsy

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MERCK

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UNIMED unimed.com.tw

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
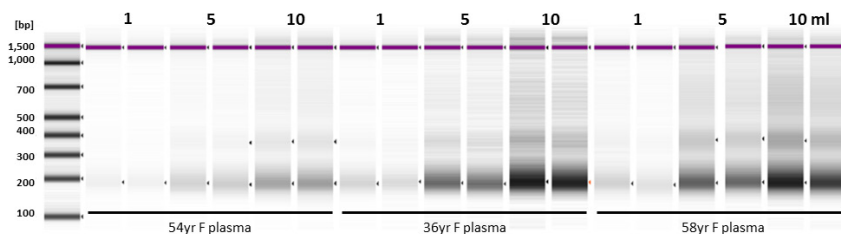
Total DNA is efficiently purified from cell-free biological fluids with the Quick-cfDNA™ Serum & Plasma Kit. Total DNA, including both high and low molecular weight species, purified (duplicates) from human maternal plasma, amniotic fluid and cerebrospinal fluid was analyzed.

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
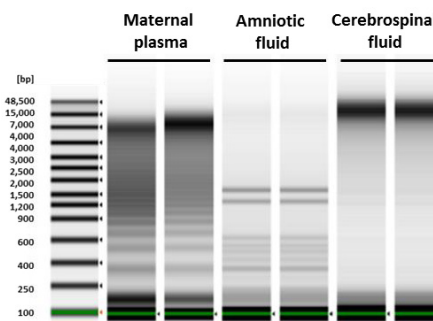
Cell-free DNA recovery scales proportionally with sample input using the Quick-cfDNA™ Serum & Plasma Kit gel image show the linear recovery of cfDNA from human plasma and serum (healthy female donors), as measured.

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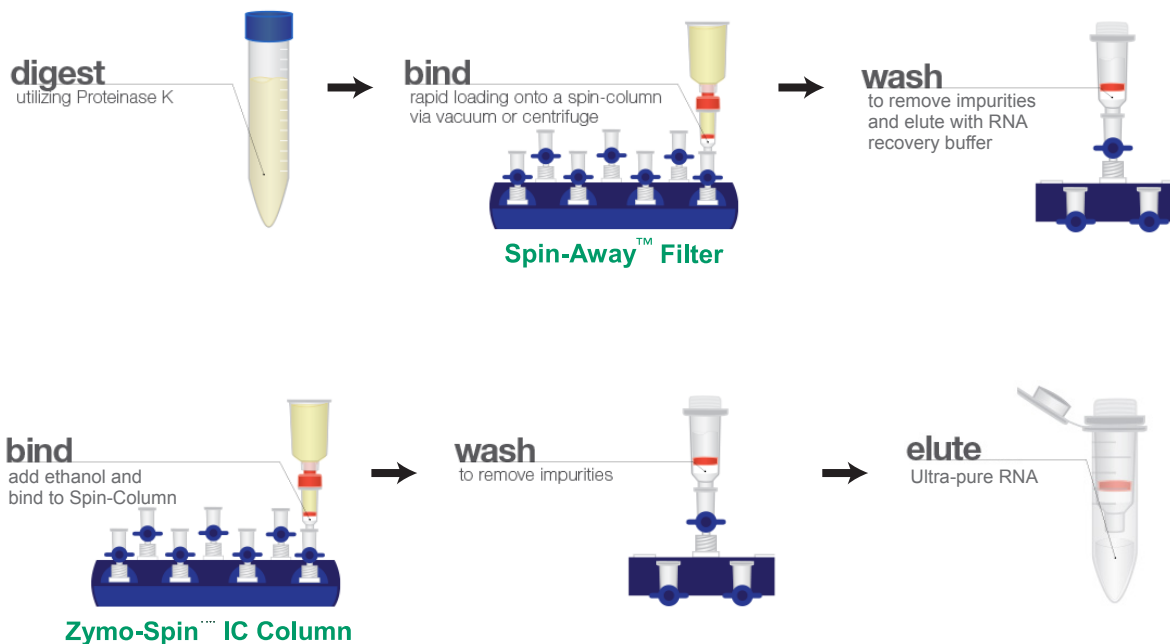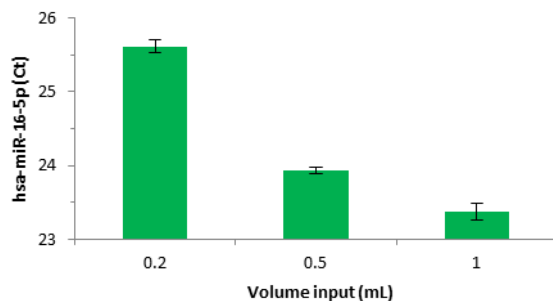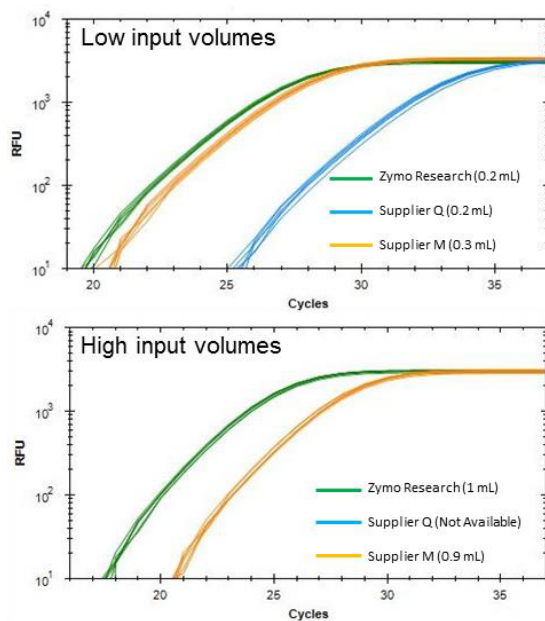
Proportional scale up relative to input volume (from 0.2 ml to 1 ml; 55y male plasma). Triplicates of purified samples used to generate RT-qPCR (Bioline; EPIK miRNA Select Kit, hsa-miR-16-5p assay).

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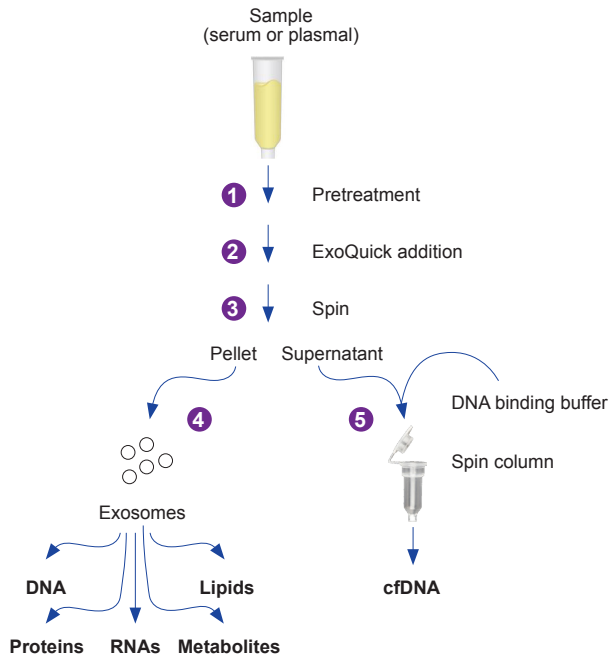
2x and 8x more yields achieved from low and high input volumes, respectively, compared to the next best product from other supplier. Triplicates of purified samples from different kits used to generate RT-qPCR data .

血漿游離 DNA (cfDNA) & 外吐小體 (Exosomes)

XCF™ COMPLETE Exosome & cfDNA Isolation Kits

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

non-exosomal protein 殘留率低

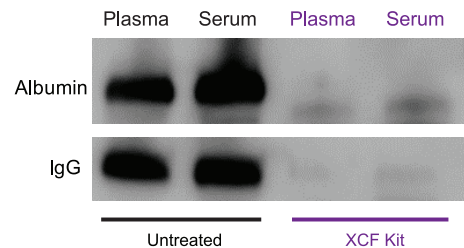The XCF COMPLETE Kit shows low protein carryover. Western blots of co-purifying human albumin (~67kD) and heavy-chain IgG (~50kD) comparing untreated serum and plasma versus the same samples processed using the XCF Kit .

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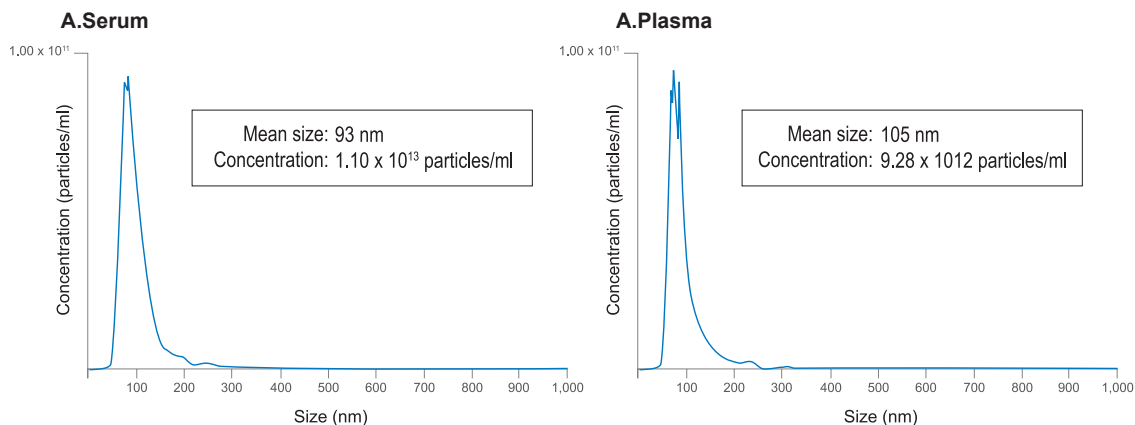
Exosomes prepared using the XCF COMPLETE Kit deliver the expected particle size profile. NanoSight NTA data for extracellular vesicles isolated from 500 μ l of serum (Panel A) and plasma (Panel B).

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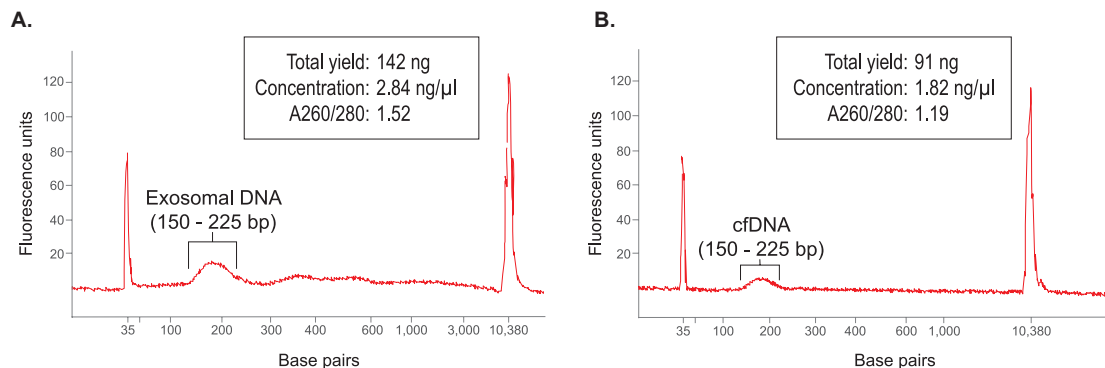搭配 **XCF™ Exosomal DNA Isolation Kit** 也可以萃取出 **exosomal DNA**

cfDNA and exosomal DNA isolated using the XCF COMPLETE Kit are of good quality and yield with the expected size. Agilent Bioanalyzer profiles of the exosomal DNA portion (Panel A) and cfDNA portion (Panel B) of 500 μ l of serum processed with the XCF COMPLETE Kit. A notable peak around ~166 bp is seen in both cfDNA and exosomal DNA.

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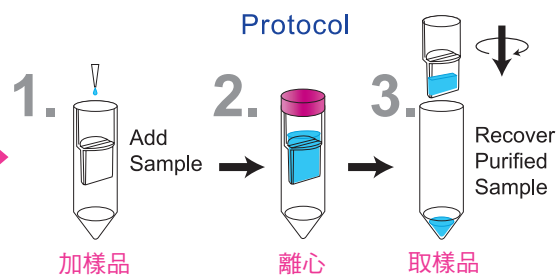
Nucleotide cut-off guidelines for single stranded DNA

Single-Stranded Nucleotide Cut-off (bp)	NMWL	
	3K (40 min)	10K (20 min)
10	98	97
15	96	95
20	94	94
25	96	96

Nucleotide cut-off guidelines for double stranded DNA

Double-Stranded Nucleotide Cut-off (bp)	NMWL		
	30K	50K	100K
137	95	95	—
301	—	94	—
657	—	99	—
1129	—	100	97

Based on >90% retention. Spinning 50 and 100K devices at higher than recommended g-force may result in lower recovery of DNA.

Recommended g-force and spin time for Amicon Ultra devices

NMWL	Maximum G-Force Rating	Spin Time (min) at 25 °C
3K	7,500	40
10K	7,500	20
30K	5,000	10
50K	2,000	20
100K	2,000	10

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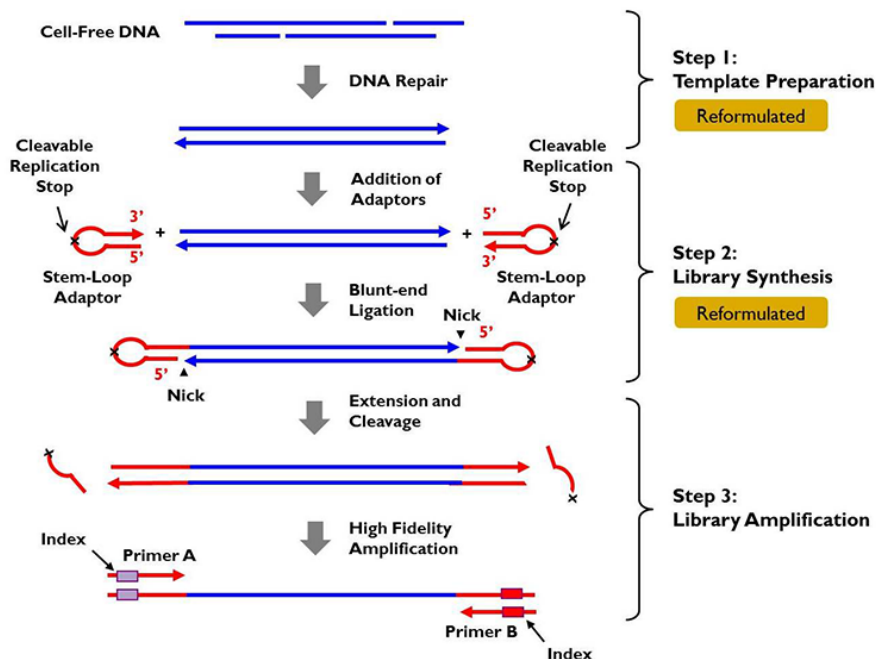
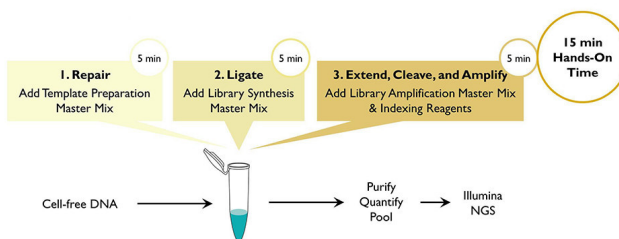
1. Cann, G.M., et al. (2012) mRNA-seq of single prostate cancer circulating tumor cells reveals recapitulation of gene expression and pathways found in prostate cancer. PLoS ONE 7(11):e49144
2. Ramsköld, D., et al. (2012) Full-length mRNA-seq from single cell levels of RNA and individual circulating tumor cells. Nature Biotechnol. 30(8):777–782.

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