MERCK

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Millipore®

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Choose the Appropriate KOD

Enzyme	KOD DNA Polymerase	KOD Hot Start DNA Polymerase	KOD XL DNA Polymerase	KOD Xtreme Hot Start DNA Polymerase
PCR Product Size (kb)	<6	<21	<30	<40
Applications	Cloning, cDNA amplification	Cloning, cDNA amplification and 70% high GC targets	Crude samples, multiplex, incorporation of derivatized dNTPs	Crude samples, long targets, difficult and 90% GC-rich targets

- KOD-XL DNA polymerase amplification results in a mixture of blunt and 3'-dA products while other KOD DNA polymerases generate blunt end products.
- NovaTaq™ is also available for routine PCR.

HOT	KOD Hot Start DNA Polymerase	71086
	KOD Hot Start Master Mix	71842
	KOD Xtreme Hot Start DNA Polymerase	71975
HOT	Agarose PCR grade	2010
	6X DNA Loading Buffer	2100
	EtBr Solution	4410
	10X TAE Buffer	574797
	10X TBE Buffer	8830
	5X TBE Buffer	8800
	Sterile Purified Water	4.86505

Bacteria Protein Exp. Cloning- pET system

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
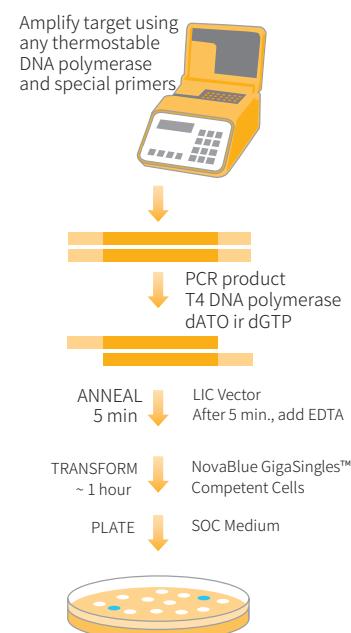
Proteinase K, lyophilized	70663 / 1.24568
Proteinase K Solution, 600 mAU/mL	71049

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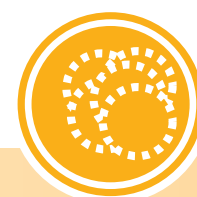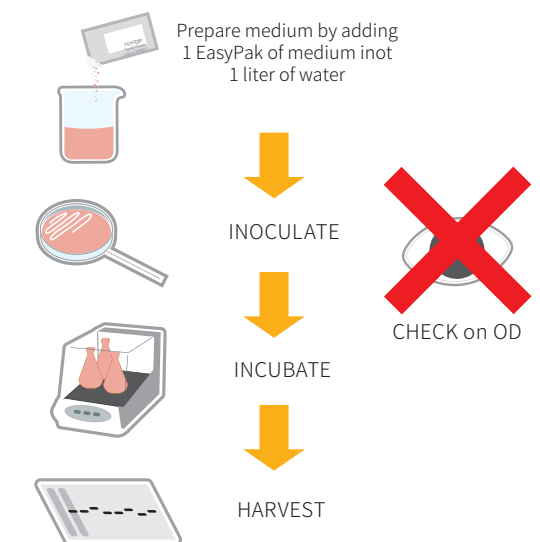
Overnight Express™ Instant LB Medium	71757-4
Overnight Express. Instant TB Medium	71491-4

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HOT	IPTG, Dioxane-Free, High Pur	420322
	X-Gal solution, 40mg/ml in DMSO	71077

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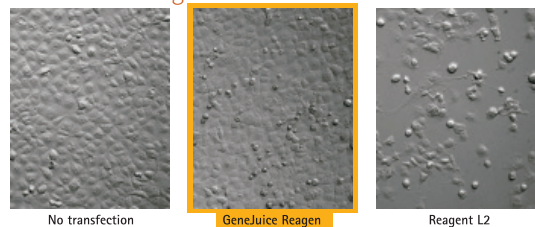
GeneJuice Reagent Is Less Cytotoxic Than Lipid-Based Transfection Reagents

GeneJuice Transfection Reagent
NanoJuice Transfection Kit
293-Free™ Transfection Reagent
ProteoJuice™ Protein Transfection
RiboJuice™ siRNA Transfection Reagent
NovaCHOice Transfection Kit

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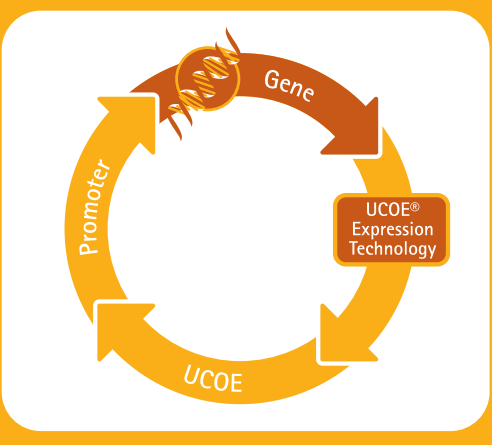
UCOE® Expression Vector, Mouse 3.2 kb, Puro 5.04865.0001
UCOE® Expression Vector, Mouse 3.2 kb, Hygro 5.04866.0001
UCOE® Expression Vector, Human, 4 kb, Puro 5.04867.0001

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When purifying proteins for functional or structural studies, the first step is to disrupt the cells or tissue sample and extract the relevant protein fraction. This step is critical because processing methods that require harsh mechanical, chemical, or enzymatic treatments can affect the target protein's integrity and activity, or otherwise expose it to degradative conditions.

Our complete range of reagents and enzymes for cell lysis and protein extraction provide you with an array of options so that you can put together the perfect extraction protocol for your particular cells and protein.

Protein Extraction Reagents Application Guide

Products by Cell Type	Starting Material		Applications			Comments
	Total Culture	Cell Pellet	1D PAGE	2D PAGE / IEF	Activity Assay	
E. coli						
BugBuster® Master Mix		■	■	■	■	Combines BugBuster® Protein Extraction Reagent with Benzonase® Nuclease and rLysozyme™ Solution. Convenient, all-in-one protein extraction reagent efficiently lyses bacteria and digests nucleic acids.
BugBuster® Protein Extraction Reagent		■	■	■	■	Efficient protein extraction from E. coli under non-denaturing conditions.
BugBuster® 10X Protein Extraction Reagent		■	■	■	■	A concentrated form of BugBuster® Protein Extraction Reagent. Ideal for extraction when a specific buffer is required for protein stability.
PopCulture® Reagent	■		■		■	Protein extraction from cells directly in the culture medium; no centrifugation required.
Yeast						
YeastBuster™ Protein Extraction Reagent		■	■		■	Efficient protein extraction from yeast under non-denaturing conditions from any volume of culture. Add 0.5 M THP Solution (included) and Benzonase® Nuclease for enhanced efficiency.
Insect						
CytoBuster™ Protein Extraction Reagent		■	■	■*	■	Gentle lysis of insect cells with retention of protein activity for assays and purification. Can use with monolayers or pellets derived from suspension cultures.
Insect PopCulture® Reagent	■		■		■	Lysis of insect cells directly in serum-free medium. Ideal for expression screening of many small samples.
Mammalian						
CytoBuster™ Protein Extraction Reagent		■	■	■*	■	Gentle lysis of mammalian cells with retention of protein activity for assays and purification. Can use with monolayers or pellets derived from suspension cultures.
ProteoExtract® Kits		■	■	■*	■	Extract protein fractions from different parts of the cell. A range of kits offering maximum flexibility.
Lysis and Extraction Enhancement						
Gram-negative bacteria (E. coli)						
rLysozyme™ Solution	■	■	■		■	Cleaves bond in peptidoglycan layer of E. coli cell wall.
Lysonase™ Bioprocessing Reagent	■	■	■		■	Convenient mixture of rLysozyme™ solution and Benzonase® Nuclease minimizes pipetting steps.
Gram-positive bacteria						
Chicken Egg White Lysozyme Solution	■	■	■		■	Cleaves bond in peptidoglycan layer of bacterial cell wall.
All cells						
Benzonase® Nuclease	■	■	■		■	Degrades all types of nucleic acids for more efficient protein extraction, faster chromatography, and reduced interference in assays.

1D PAGE — One-dimensional Polyacrylamide Gel Electrophoresis

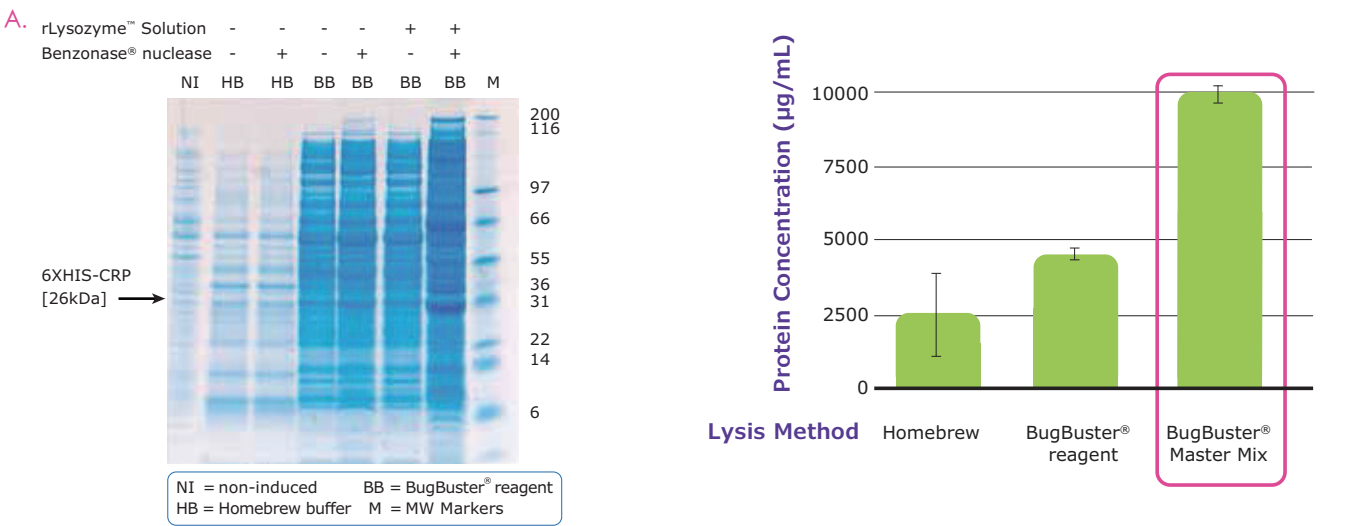
2D PAGE — Two-dimensional Polyacrylamide Gel Electrophoresis

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BugBuster® Protein Extraction Kits and Reagents

BugBuster® reagent is superior to “homebrew” lysis buffer and BugBuster® reagent with both Benzonase® nuclease and rLysozyme™ solution produced lysates with the highest 6XHIS-CRP yields. (A) E. coli lysates (5 µL of 1 mL total lysate) from various lysis protocols were fractionated and analyzed by SDS-PAGE. A band

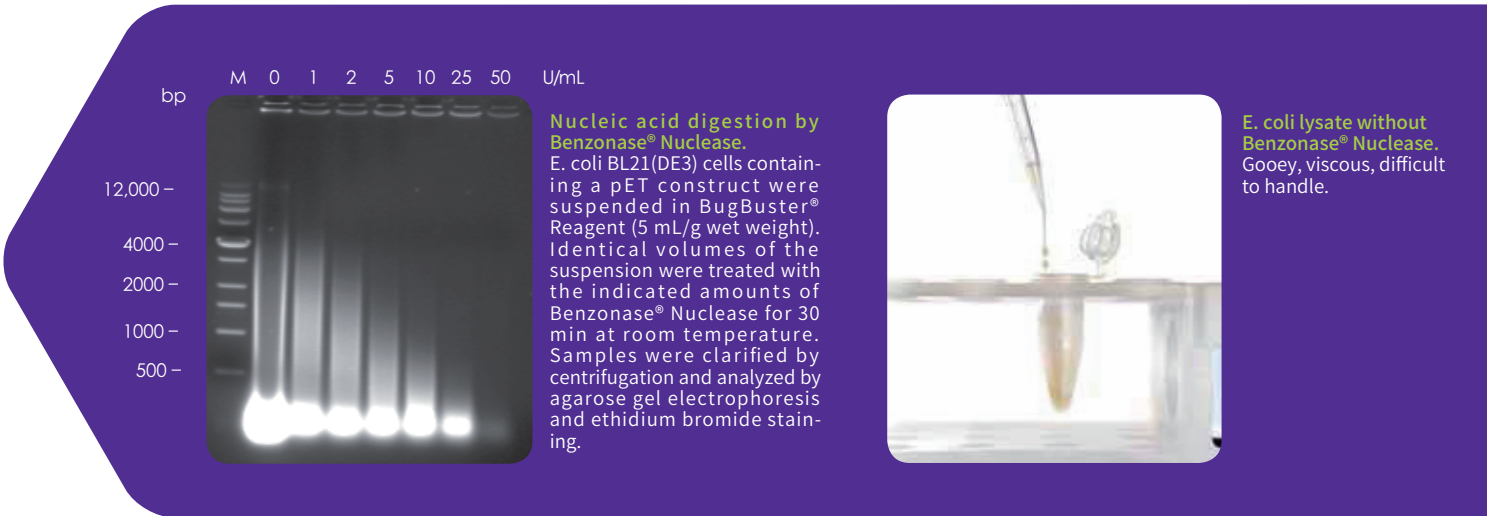
corresponding to 6XHIS-CRP is prominently visualized in the BB +/- lane. (B) Cleared cell lysates (2 µL of 1 mL total) were spotted on assay cards and quantified using the Direct Detect® spectrometer. In each case, bars represent the average of 3 independent samples.

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Benzonase® Nuclease
Effectively reduce viscosity and remove nucleic acids from protein solutions
Benzonase® Nuclease is a genetically engineered endonuclease from *Serratia marcescens*. It degrades all forms of DNA and RNA (single stranded, double stranded, linear and circular) while having no proteolytic activity. It is effective over a wide range of conditions and has an exceptionally high specific activity. Benzonase® nuclease is an excellent choice for viscosity reduction to shorten processing time and increase protein yields.

- Benzonase® Advantages**
- Compliant with FDA guidelines for nucleic acid contamination
 - Functional between pH 6 and 10, from 0 °C to 42 °C, for maximum versatility
 - Active in the presence of ionic and non-ionic detergents, reducing agents, PMSF (1 mM), EDTA (1 mM) and urea.
 - Available in ultrapure (> 99 % by SDS-PAGE) and pure (> 90 %) grades
 - Available in standard concentration (25 U/µL) and high concentration (HC, 250 U/µL).

CHEMICON®’s Nuclear Extraction Kit (Catalog No. 2900) provides a simple and convenient method for the isolation of cytoplasmic and nuclear samples from mammalian cell culture or tissue samples. The Nuclear Extraction Kit can be used in the preparation of purified proteins for use in Western blotting, Electrophoretic Mobility Shift Assays (EMSA), and in CHEMICON®’s Transcription Factor Assay product line.


Ordering Information

Application	Description	Catalog No.
Bacteria	BugBuster® Protein Extraction Reagent	70584
	BugBuster® Master Mix	71456
	BugBuster® Plus Benzonase® Nuclease	70750
	BugBuster® Plus Lysonase™ Kit	71370
	BugBuster® 10X Protein Extraction Reagent	70921
	PopCulture® Reagent	71092
Mammalian	CytoBuster™ Protein Extraction Reagent	71009
	NucBuster™ Protein Extraction Reagent	71183
	PhosphoSafe™ Extraction Reagent	71296
Yeast	YeastBuster™ Protein Extraction Reagent	71186
Insect	Insect PopCulture® Reagent	71187

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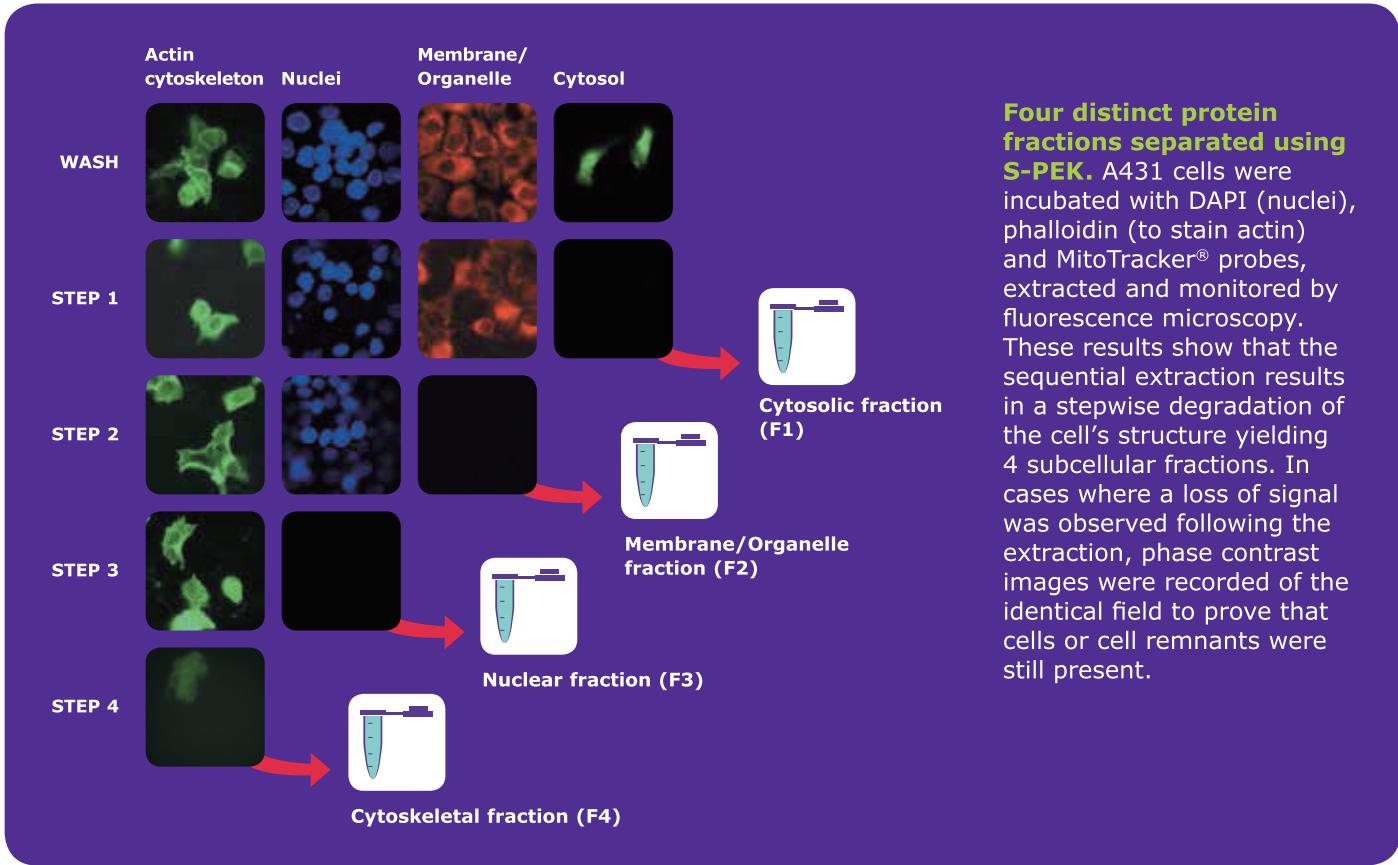
ProteoExtract® Subcellular Proteome Extraction Kit (S-PEK)
Reproducible extraction of subcellular proteomes from mammalian cells.
539790 - IKITCN

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ProteoExtract® Native Membrane Protein Extraction Kit (M-PEK)
Isolation of native membrane proteins from mammalian cells and tissue.
Extract proteins associated with cellular membranes with M-PEK. Extremely mild extraction conditions yield a 3–5 fold enrichment of integral membrane and membrane-associated proteins. The simple, two-step procedure enables processing of multiple samples in parallel. Extraction from tissues requires isolation of viable cells before proceeding with the extraction protocol.

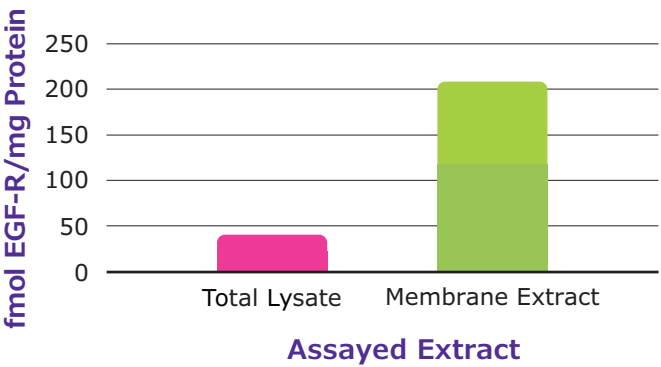Applications for Extracted Membrane Proteins:

Notably increased enrichment of EGF receptor using M-PEK compared to total cell lysate. HEK293 cells were extracted with buffered 1 % Triton® X-100 surfactant to generate a total lysate or extracted with M-PEK to yield a membrane fraction. Equal volumes of these fractions were utilized to quantitate the concentration of EGF receptor in the samples using an EGF-R ELISA Kit. Protein concentrations were used to calculate the amount of EGF-R per mg protein in the total lysate and the membrane fraction. The measurements demonstrate a 4.5 fold enrichment of the EGF receptor in the M-PEK-extracted membrane fraction.

Ordering Information

Application	Description	Catalog No.
Organelle Fractionation	ProteoExtract® Subcellular Protein Extraction Kit	539790
	ProteoExtract® Complete Mammalian Protein Extraction Kit	539779
	ProteoExtract® Cytosol/Mitochondria Fractionation Kit	QIA88
	ProteoExtract® Native Cytoskeleton Enrichment Kit	17-10210
	ProteoExtract® Cytoskeleton Enrichment and Isolation Kit	17-10195
Membrane Proteins	ProteoExtract® Native Membrane Protein Extraction Kit	444810
	ProteoExtract® Transmembrane Protein Extraction Kit	71772
Mass Spec Peptide Enrichment	ProteoExtract® All-in-One Trypsin Digestion Kit	650212
	ProteoExtract® Glycopeptide Enrichment Kit	72103
	ProteoExtract® Phosphopeptide Enrichment TiO ₂ Kit	539722

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買任何 Merck 產品
皆可以加價購 1ml
Protease Inhibitor Cocktail III
(#539134, ready-to-use, 100X)

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買任何 Merck 產品
皆可以加價購 1ml
Phosphatase Inhibitor Cocktail V
(#524629, ready-to-use, 50X)

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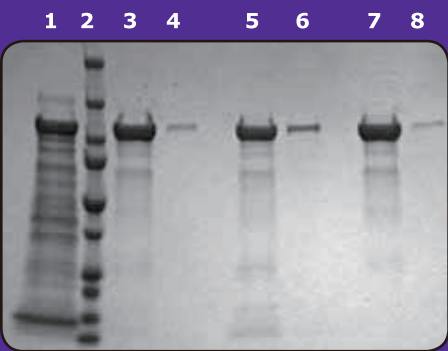
Protease Inhibitor Cocktail III +
Phosphatase Inhibitor Cocktail V
各1ml 包裝
讓您的蛋白萬無一失！

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His•Tag® Purification

Ni-NTA His•Bind® Resin has a binding capacity of over 10 mg of His-Tagged fusion protein per mL resin. The agarose matrix on the Ni-NTA His•Bind® Superflow™ Resin is structured with more crosslinking for enhanced bead rigidity, for exceptional compatibility with FPLC. Our IDA His•Bind® resins are offered uncharged to allow flexibility of choice in the metal ion (Nickel, Cobalt, Zinc, Iron, Copper, etc.). IDA supports can be recycled many times with no loss in performance.

Lane	Sample
1	Crude Extract
2	Markers
3	Ni-NTA Competitor Q Elution
4	Ni-NTA Competitor Q Strip
5	Ni-NTA Competitor G Elution
6	Ni-NTA Competitor G Strip
7	Ni-NTA His•Bind® Elution
8	Ni-NTA His•Bind® Strip


Ni-NTA His•Bind® performance vs. equivalent competitor resins Vector pET-28b (+) was used to express a His-Tag fusion protein of 119KDa in *E. coli* BL21 (DE3) cells, induced culture was processed with BugBuster® Master Mix, and protein extract was divided evenly to proceed to the His-Tag purification using Ni-NTA His•Bind®, Ni-NTA Competitor Q and Ni-NTA Competitor G resins. Ni-NTA His•Bind® resins show higher binding capacity and a better purification.

Ordering Information

Application	Description	Catalog No.
Ni-NTA His•Bind® Resin		
Small to medium scale	Ni-NTA His•Bind® Resin	70666
Gravity flow column	BugBuster® Ni-NTA His•Bind® Purification Kit	70751
Recommended for eukaryotic extracts	Ni-NTA Buffer Kit	70899
Ni-NTA His•Bind® Superflow™ Resin		
Small to production scale	Ni-NTA His•Bind® Superflow™ Resin	70691
FPLC or gravity flow column	Ni-NTA Buffer Kit	70899
Uncharged IDA His•Bind® Resin		
Uncharged (metal flexibility)	IDA His•Bind® Resin	69670
Reusability	His•Bind® Buffer Kit	69755
Small to medium scale	His•Bind® Purification Kit	70239
Gravity flow column or batch mode	BugBuster® His•Bind® Purification Kit	70793

GST•Tag™ Purification

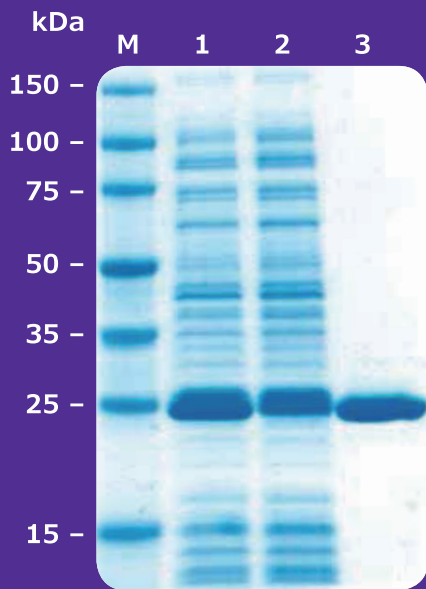The GST fusion system is based on the widely recognized affinity of glutathione-S-transferase (GST) fusion proteins for immobilized glutathione. Our GST Resin utilizes an 11-atom spacer arm to covalently attach reduced glutathione to the solid support via a sulfide linkage. The resin can be reused several times without loss of capacity, and the high degree of substitution of glutathione ensures exceptional binding capacity.


Lane	Sample
M	PerfectProtein™ markers 15–150 kDa
1	BugBuster® extract
2	Flow-through
3	Eluate

GST•TAG™ Purification
GST•Bind™ purification. A crude extract containing unfused GST was applied to a 2 mL GST•Bind™ Resin column. Total protein yield after purification was 8 mg/mL resin.

S•Tag™ Purification

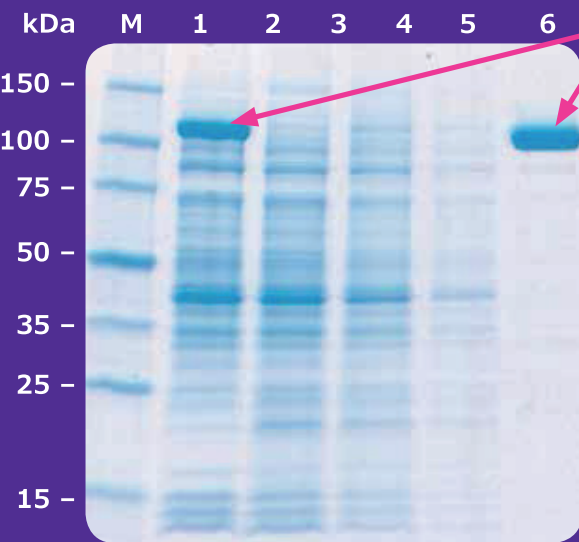The S•Tag™ fusion protein is a short 15-aa sequence that specifically binds with high affinity the 104-aa S-Protein ($K_D = 10^{-9}$ M, 1000 times stronger than the interaction between nickel and His•Tag® fusion protein). Fusion proteins can be easily purified by cleavage with site-specific proteases or in acidic buffers.

Lane	Sample
M	PerfectProtein™ markers 15–150 kDa
1	Crude extract
2	Flow-through
3	Wash 1
4	Wash 2
5	Eluate + Biotinylated Thrombin
6	Eluate after Biotinylated Thrombin removal

S•Tag™ affinity purification
S•Tag™ β-gal expressed from a pET construct was purified from a crude soluble fraction using S-protein Agarose under native conditions. Elution of the target protein from the agarose was performed by digestion with Biotinylated Thrombin, which was subsequently removed with Streptavidin Agarose. The fractions are indicated.

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Restriction and Biotinylated Grade **Thrombin**

Highly efficient, specific cleavage of fusion proteins

Restriction Grade Thrombin is qualified to specifically cleave target proteins containing the recognition sequence LeuValProArgâGlySer. The preparation is functionally tested for activity with fusion proteins and is free of detectable contaminating proteases. Thrombin is supplied with 10X Thrombin Cleavage Buffer and a Cleavage Control Protein. Biotinylated Thrombin is identical in activity to Restriction Grade Thrombin, but has covalently attached biotin for easy removal of the enzyme from cleavage reactions using immobilized streptavidin. Our Thrombin Cleavage Capture Kit includes not only biotinylated thrombin and immobilized streptavidin, but also all required buffers and filters for complete, convenient recovery of cleaved protein.

HRV 3C Protease

Highly efficient, specific cleavage of fusion proteins

Recombinant type 14 3C protease from human rhinovirus (HRV 3C) is a highly purified, recombinant 6XHis-tagged enzyme, which recognizes the cleavage site LeuGluValLeuPheGlnâGlyPro. The small, 22 kDa size of the protease, with optimal activity at 4 °C, high specificity, and His-tag fusion make HRV 3C protease an ideal choice for rapid removal of fusion tags.

Strep · Tag® II Purification

The Strep · Tag® fusion protein II is an 8 amino acid sequence that binds to the biotin pocket of Streptavidin with 100 times higher binding capacity.

T7 · Tag® Purification

Purification is antibody-based. Covalently coupled to agarose beads, the T7 · Tag® monoclonal antibody captures the T7 · Tag® epitope — a sequence of 11

Streptavidin Agarose

Cross-linked agarose is covalently coupled with pure streptavidin under controlled conditions. The stable linkage to the resin minimizes leaching of the streptavidin while maintaining full binding activity.

Ordering Information

Description	Catalog No.
Restriction-Grade Thrombin	69671
Biotinylated Thrombin	69672
Thrombin Cleavage Capture Kit	69022
Restriction Grade Factor Xa	69036
Factor Xa Cleavage Capture Kit	69037
Recombinant Enterokinase	69066
Enterokinase Cleavage Capture Kit	69067
HRV 3C Protease	71493
Tag•off™ High Activity rEK	71537
Tag•off™ rEK Cleavage Capture Kit	71540

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
High speed immunoprecipitation with magnetic beads compared to agarose. In parallel indirect immunoprecipitations, PureProteome™ magnetic beads offered a 50 % reduction in incubation time while yielding results equivalent to agarose beads.

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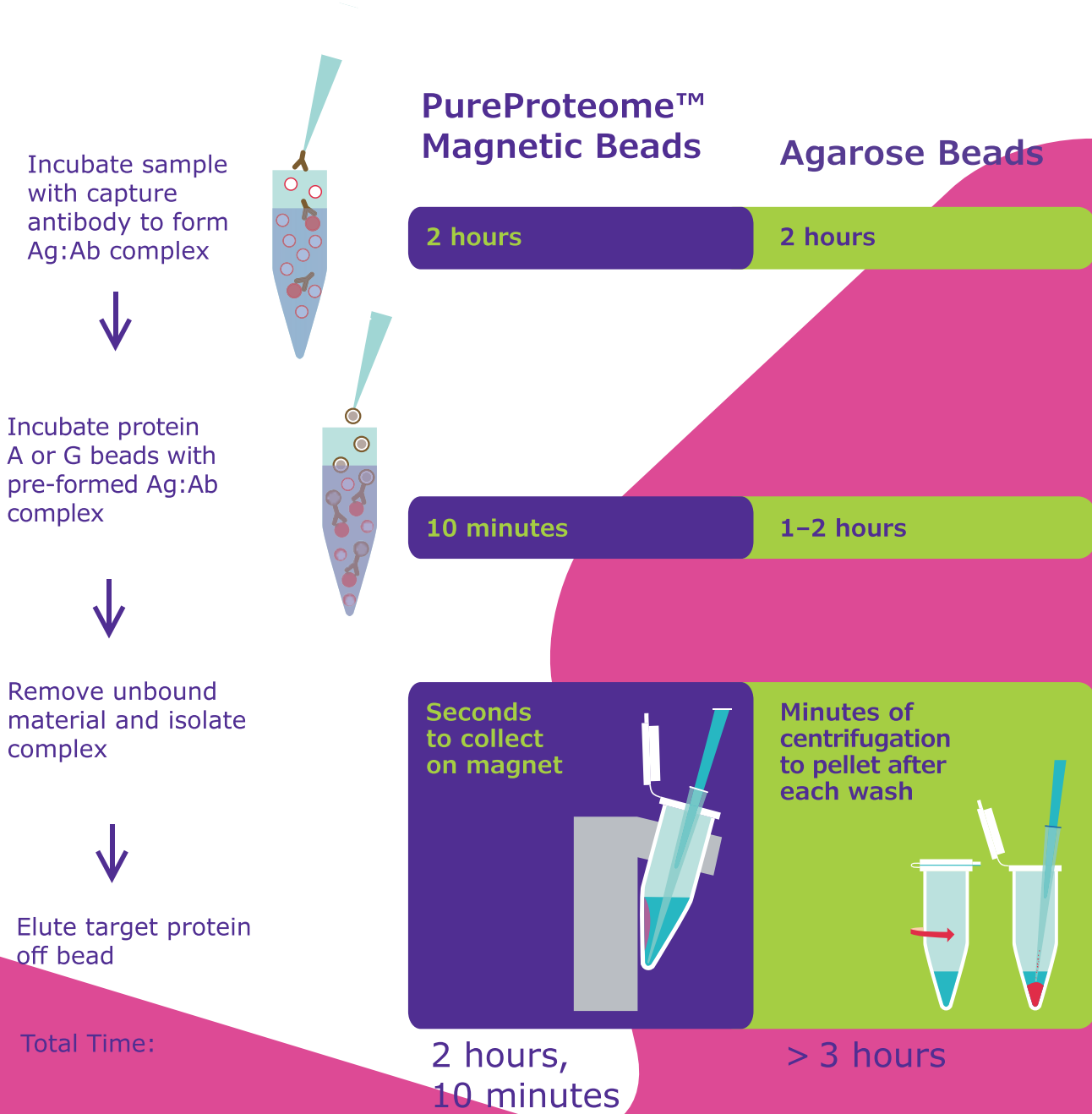Ordering Information

Application	Description	Catalog No.
IP, Antibody Purification, F(ab) Purification	PureProteome™ Protein A Magnetic Beads	LSKMAGA10
	PureProteome™ Protein G Magnetic Beads	LSKMAGG10
	PureProteome™ Protein A/G Mix Magnetic Beads	LSKMAGAG10
	PureProteome™ Kappa Ig-Binder Magnetic Beads*	LSKMAGKP02
	PureProteome™ Lambda Ig-Binder Magnetic Beads*	LSKMAGLM02
Biotinylated Molecule Purification	PureProteome™ Streptavidin Magnetic Beads	LSKMAGT10
His•Tag® Tagged Protein Purification	PureProteome™ Nickel Magnetic Beads	LSKMAGH10
Custom Labelled (Flexibility to Bind Ligand of Choice)	PureProteome™ NHS FlexiBind Magnetic Beads	LSKMAGN04
	PureProteome™ Carboxy FlexiBind Magnetic Beads**	LSKMAG1CBX10
Magnetic Stands	PureProteome™ Magnetic Stand, 8-well	LSKMAGS08
	PureProteome™ Magnetic Stand, 15 mL	LSKMAGS15

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IP-05 BSA blocking proteinA/G agarose beads

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快試試 BSA block beads!

Magna Grip™ Rack contains 8 holes suitable for 1.5-2.0 ml tubes or spin columns. May also be used with either 15 ml or alternately, 0.5 ml tubes, simply by standing the rack on either side.

Designed for immunoprecipitation applications. This product is blocked with BSA to reduce non-specific binding and cannot be used for purification.

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ProteoExtract® Agarose Columns

Human serum and plasma samples are rich sources of proteomic information, reflecting processes regulating normal or diseased states. Today’s ultra-sensitive analytical methods, such as two dimensional (2D) gel electrophoresis and mass spectrometry, can detect minute changes in expression profiles — but ultrasensitive approaches typically require the removal of highly abundant proteins (HAP) and moderately abundant proteins (MAP).

We offer a range of kits and resins for depleting high-abundance proteins (HAP) from serum or plasma samples. Choose from the PureProteome™ magnetic bead kits and resins or the ProteoExtract® agarose columns. First, identify the species of your serum/plasma source — the following tables summarize the different solutions for your needs.

Ordering Information

Description	Format	Species	Proteins Depleted	Catalog No.
ProteoExtract® Albumin/IgG Depletion Kit	Agarose columns	Human, rabbit, rat, mouse, pig and bovine	Albumin and IgG > 80 %	122642
ProteoExtract® Albumin Depletion Kit			Albumin > 80 %	122640

Application	Description	Catalog No.
Depletion/Enrichment	PureProteome™ Albumin Magnetic Beads	LSKMAGL10
	PureProteome™ Albumin/IgG Depletion Kit	LSKMAGD12
	PureProteome™ Human Albumin/Immunoglobulin Depletion Kit*	LSKMAGHDKIT

*Human only.

ProteoPrep® Immunoaffinity Albumin & IgG Depletion Kit

The ProteoPrep® Immunoaffinity medium in pre-packed spin columns is a mixture of two-beaded media containing recombinantly expressed, small single-chain antibody ligands, providing low non-specific binding and high capacity. This kit is also effective in depleting albumin and IgG from mouse and guinea pig serum.

The ProteoPrep® Blue Albumin & IgG Depletion Kit is designed to specifically remove albumin and IgG from 25 samples of human serum (25 µL to 100 µL) in preparation for two-dimensional electrophoresis (2DE). Albumin-depleted serum samples generated by this kit are in urea rather than salt-based buffers with little dilution, eliminating the need to precipitate the sample prior to performing 2DE.

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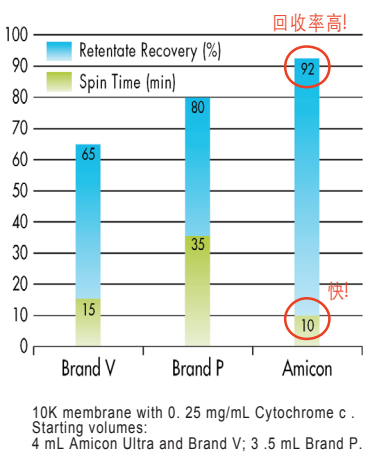
raw material 濃縮與 Protein Fractionation, 不用擔心樣品流失

Amicon Ultra-4 Centrifugal Filter

Protein (Concentration)	NMWL*	Retentate Recovery (%) by Brand			
		V 6 mL	PG	P 7 mL	Amicon Ultra-4
Cytochrome C (0.025 mg/mL)	5,000	13	—	—	96
	9,000	—	—	56	—
	10,000	0	9	—	97
BSA (0.1 mg/mL)	20,000	—	—	44	—
	30,000	76	89	—	96
	50,000	0	—	—	79
IgG (0.1 mg/mL)	100,000	78	76	—	96

Protein (Concentration)	NMWL	Dilute Protein Recovery (%) by Brand			
		V 6 mL	P 7 mL	Amicon Ultra-4	
BSA (100 µg/mL)	30,000	76	80	96	
BSA (10 µg/mL)	30,000	51	66	93	
BSA (1 µg/mL)	30,000	16	67	88	
BSA (0.14 µg/mL)	30,000	6	44	87	

*Nominal Molecular Weight Limit or membrane cut-off

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1. What are you concentrating?
2. What is the source of your sample?
3. What is the molecular weight of your protein?
4. What is your starting volume?
5. What is your desired final volume?

Amicon Ultra Selector Tool

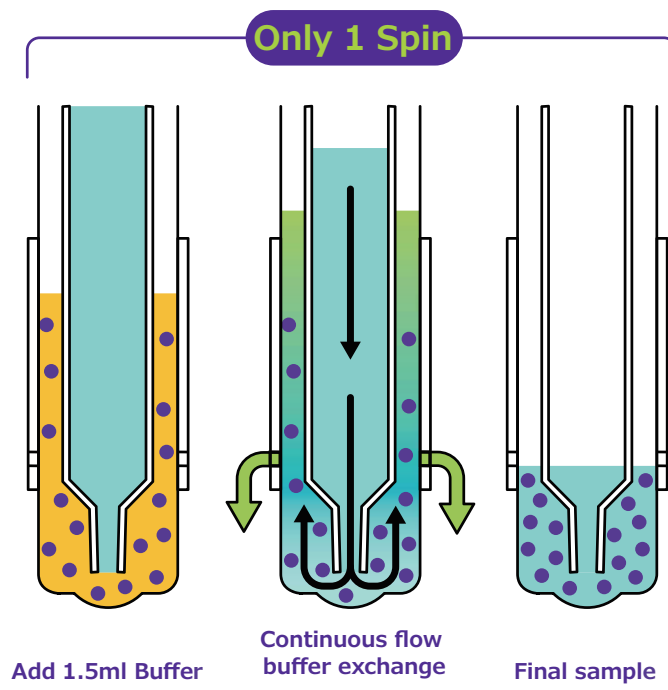
FAST AND EASY CONCENTRATION

What are you concentrating?

- ☐ DNA
- ☒ Protein or other molecules of known weight in kilodaltons (kDa)
- ☐ Mammalian cells
- ☐ Bacterial cells
- ☐ Virus
- ☐ Other molecules/particles of known size in nanometers

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Ordering Information

To choose the appropriate Amicon® Pro device, determine the nominal molecular weight cut-off (NMWCO) for your protein of interest and your desired affinity purification scheme.

Amicon® Pro purification system — No Reagents Included	NMWCO				
	3,000	10,000	30,000	50,000	100,000
Amicon® Pro Purification System 12/pk	ACS500312	ACS501012	ACS503012	ACS505012	ACS510012
Amicon® Pro Purification System 24/pk	ACS500324	ACS501024	ACS503024	ACS505024	ACS510024
Amicon® Pro Purification System 24/pk without Amicon® Ultra 0.5 mL filter			ACS500024		

大體積濃縮交給 AMicon stirred cell

Amicon® Stirred Cells

50 mL to 400 mL concentration

Amicon® stirred cells provide high flow rates with solutions up to 10 % macrosolute concentration and are capable of rapid concentration, or salt removal followed by concentration in the same unit. For protein concentration, gas pressure is applied directly to ultrafiltration cell. Solutes above the membrane’s nominal molecular weight cut-off (NMWCO) are retained in cell, while water and solutes below the cut-off pass into the filtrate and out of cell.


Advantages

- Gentle magnetic stirring minimizes concentration polarization and shear denaturation.
- All stirred cells can be autoclaved.
- Three different sizes to handle volumes from 50 mL to 400 mL
- High flow rates with solutions up to 10 % macrosolute concentration

Available in three sizes

Max. Working Volume	Catalog No.
50 mL	UFSC05001
200 mL	UFSC20001
400 mL	UFSC40001

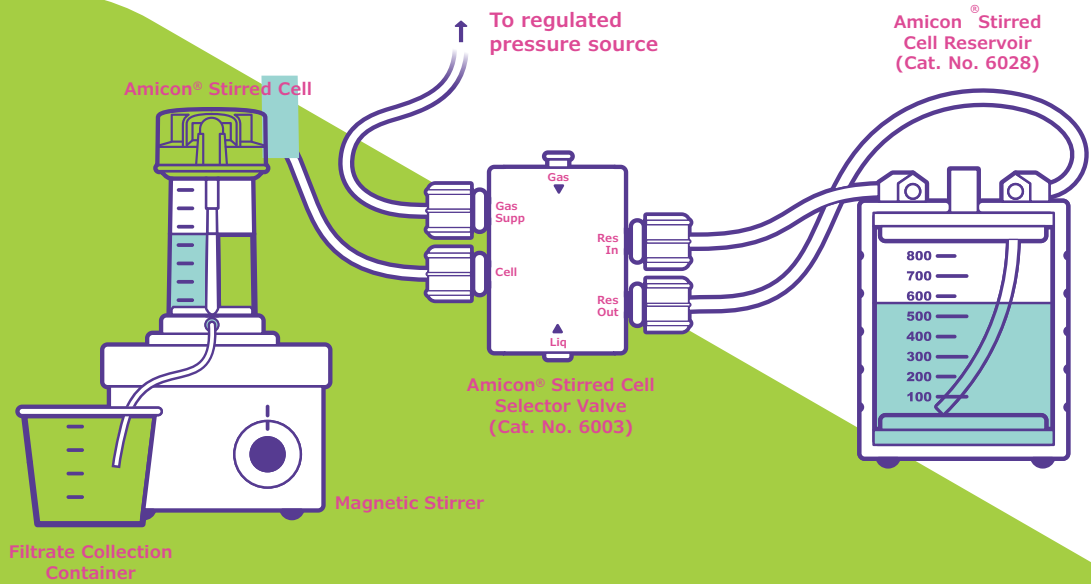
Amicon® Stirred Cell setup for large volume concentration and continuous diafiltration using selector valve and reservoir accessories.

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	Description	Cat. No.	Size
Organics & Inorganics	OmniPur® Agarose	2120	100 g
	OmniPur® Agarose	2125	500 g
	OmniPur® Agarose, Low Melting	2070	100 g
	OmniPur® Glycerol	4750	1 l
	OmniPur® Glycerol	4760	4 l
	OmniPur® Phenylmethyl Sulfonyl Fluoride	7110	5 g
	OmniPur® Phenylmethyl Sulfonyl Fluoride	7130	250 g
	OmniPur® Protein Molecular Weight Marker	7331	500 uL
	OmniPur® TEMED	8920	25 ml
Buffers	OmniPur® Glycine	4810	500 g
	OmniPur® Glycine	4840	5 kg
	OmniPur® TRIS	9210	500 g
	OmniPur® TRIS	9230	5 kg
	OmniPur® TRIS Hydrochloride	9310	500 g
	OmniPur® TRIS Hydrochloride	9350	5 kg
Detergents	OmniPur® Sodium Dodecyl Sulfate, SDS	7910	500 g
	OmniPur® Tween® 20 (Polyoxyethylene (20) Monolaurate)	9480	1 l
	OmniPur® Tween® 20 (Polyoxyethylene (20) Monolaurate)	9485	4 l
	OmniPur® Tween® 80 (Polyoxyethylene (20) Monooleate)	9490	1 l
	OmniPur® Tween® 80 (Polyoxyethylene (20) Monooleate)	9495	4 l
	OmniPur® Triton® X-100 Surfactant	9400	100 ml
	OmniPur® Triton® X-100 Surfactant	9440	4 l
	OmniPur® Triton® X-100 Surfactant	9460	20 l
Salts, Acids & Bases	OmniPur® Sodium Chloride	7710	500 g
	OmniPur® Sodium Chloride	7760	5 kg
	OmniPur® Magnesium Chloride, Hexahydrate	5980	500 g
	OmniPur® Ammonium Persulfate	2300	25 g
Buffers	OmniPur® 10X PBS Liquid Concentrate	6505	4 l
	OmniPur® 10X PBS Liquid Concentrate	6504	20 l
	OmniPur® 10X PBS Liquid Concentrate	6507	4 l
	OmniPur® 10X PBS, Premixed Powder	6508	1 l
	OmniPur® PBS Tablets	6501	200 ea
	OmniPur® 10X TBE Buffer	8820	4 l
	OmniPur® 10X TBE Buffer	8830	20 l
	OmniPur® 10X TBE Buffer	8840	200 l
	OmniPur® 10X TG Buffer	9000	4 l
	OmniPur® 10X TG-SDS Buffer	9015	4 l
Chelating Agents	OmniPur® EDTA	4005	500 g
	OmniPur® EDTA	4008	5 kg
	OmniPur® EDTA, Disodium Salt, Dihydrate	4010	250 g
Dyes and Stains	OmniPur® 6X Loading Buffer	2100	5ml
	OmniPur® Bromophenol Blue, Sodium Salt	2830	25 g
	OmniPur® Ethidium Bromide	4410	10ml
Water	Water, WFI Quality, Sterile Purified Water	4.86505.0500	500 ml
	Water, WFI Quality, Sterile Purified Water	4.86505.1000	1 l
	Water, WFI Quality, Sterile Purified Water	4.86505.5000	5 l

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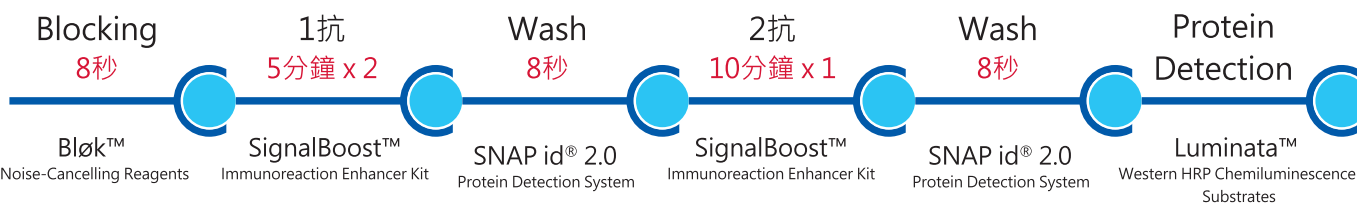reblot™ plus western blot recycling

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SNAP i.d.® Basic Package

Components	Package Catalogue No.
SNAP i.d.® 2.0 Protein Detection System – please specify which size	BASICMINI
Immobilon®-P Membrane, 26.5 cm x 3.75 m roll	BASICMIDI
SNAP i.d.® 2.0 Blot Holders, 100 units/pk	BASICMM

Western Blotting Package

Components	Package Catalogue No.
Immobilon®-P Membrane, 26.5 cm x 3.75 m roll	WBPRO
Bløk® Noise Cancelling Reagent	
Luminata™ Classico Western HRP Substrate, 100 mL	
Reblot™ Plus Strong Antibody Stripping Solution, 10x, 50 mL	

Fluorescent Blotting Package

Components	Package Catalogue No.
Immobilon®-FL PVDF Membrane, 0.45 µm, 26.5 cm x 3.75 m roll	WBPROFL
Bløk® Noise Cancelling Reagent	

Luminata™ Starter Package

Components	Package Catalogue No.
Luminata™ Classico Western HRP Substrate, 100 mL	WBLUM0100
Luminata™ Crescendo Western HRP Substrate, 100 mL	
Luminata™ Forte Western HRP Substrate, 100 mL	

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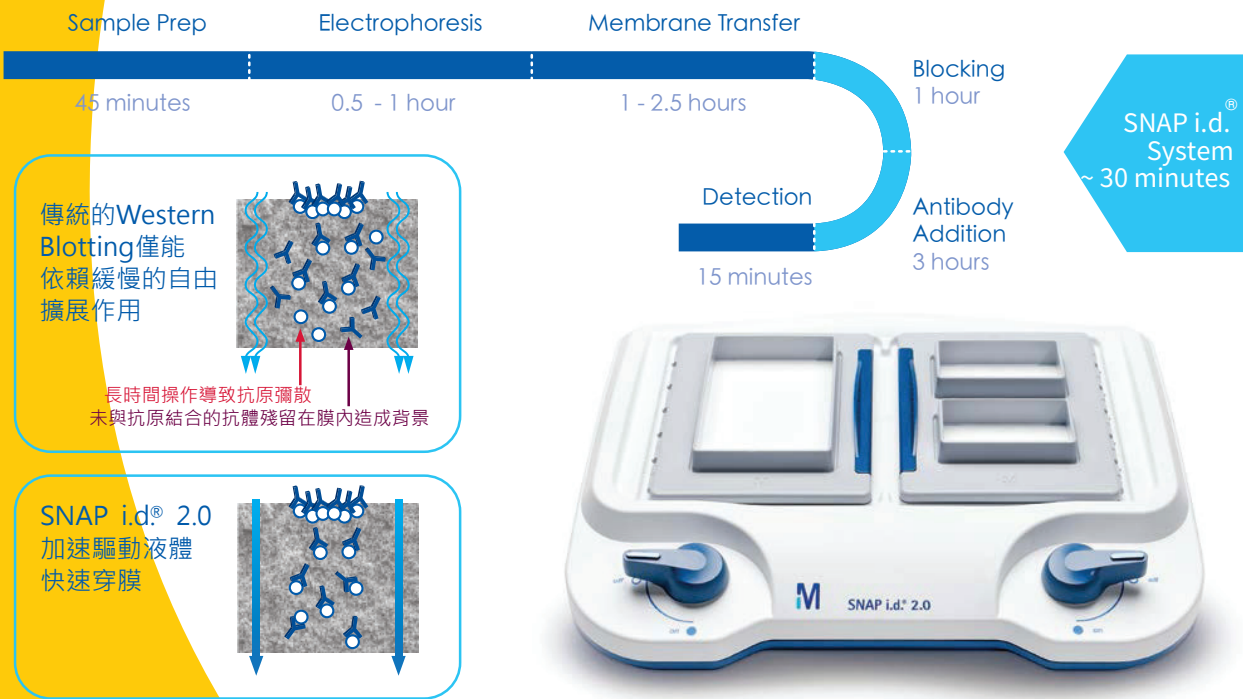
SNAP i.d.® 2.0
 Protein Detection System for Western Blotting

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	Luminata™ Classico	Luminata™ Crescendo	Luminata™ Forte
Unique Feature	Premixed	Premixed	Premixed
Best used for	Blots where the primary antibody is incubated for ~1 hr	Blot where the primary antibody is incubated > 2 hrs	Blots with overnight primary antibody incubation, or detection of PTM** proteins
Detection Range	~6 pg	~1–3 pg	~400 fg
Signal Duration	1 hr	3 hr	3 hr
Stock Solution Stability	1 yr at 4 °C	1 yr at 4 °C	1 yr at room temperature

**PTM - Post-translationally modified.

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
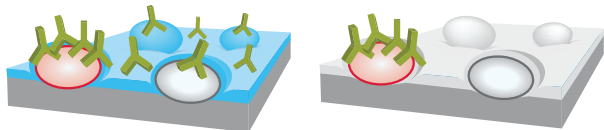
IMMOBILON PVDF

How do Immobilon® membranes work?

Membranes bind biomolecules through hydrophobic (polyvinylidene fluoride (PVDF) or electrostatic (cellulose-based membranes) interactions. Membrane pores increase the surface binding area while restricting sizes of bound proteins.

Key Benefits

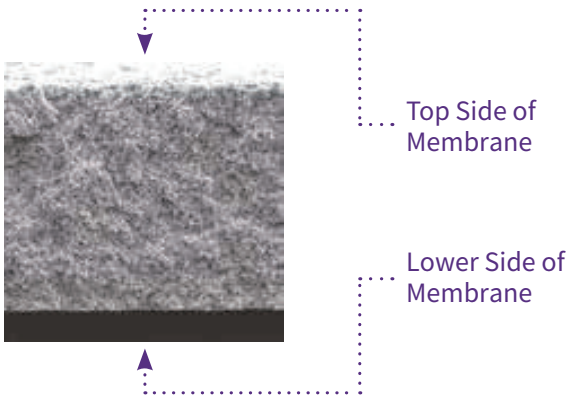- Stronger protein signals due to high adsorption & retention of protein in the membrane
- Prolonged shelf life due to higher tensile strength
- Easier stripping & reprobing with PVDF membranes
- A variety of pore sizes provides optimal protein retention

Membranes are 3-dimensional structures full of microscopic pores (Scanning electron microscope image of a crosssection of Immobilon®-P Membrane, Magnification: 500x).

Membrane comparison table

	PVDF Membranes				Nitrocellulose Membranes			
Manufacturer	Millipore® Brand				GE Healthcare Life Sciences			
Membrane Name	Immobilon®-E	Immobilon®-P	Immobilon®-FL	Immobilon®-P ⁹⁰	Amersham® Protran	Amersham® Protran premium	Amersham® Protran	Amersham® Protran premium
Chemistry	PVDF	PVDF	PVDF	PVDF	Nitrocellulose	Nitrocellulose	Nitrocellulose	Nitrocellulose
Pore Size	0.45 µm	0.45 µm	0.45 µm	0.2 µm	0.45 µm	0.45 µm	0.2 µm	0.2 µm
Detection methods	Chemiluminescence, Colorimetric, Fluorescence	Chemiluminescence, Colorimetric	Chemiluminescence, Colorimetric, Fluorescence (including near-infrared)	Chemiluminescence, Colorimetric, Fluorescence	Chemiluminescence, Colorimetric	Chemiluminescence, Colorimetric, Fluorescence	Chemiluminescence, Colorimetric	Chemiluminescence, Colorimetric, Fluorescence
Goat IgG	225-293 µg/cm²	294 µg/cm²	300 µg/cm²	448 µg/cm²	115-125 µg/cm²	162-180 µg/cm²	150-176 µg/cm²	173-203 µg/cm²
BSA	N/A	215 µg/cm²	205 µg/cm²	340 µg/cm²	N/A	N/A	N/A	N/A
Insulin	N/A	160 µg/cm²	155 µg/cm²	262 µg/cm²	N/A	N/A	N/A	N/A
Applications	Western blots	Western blots, dot blots			Western, Southern and Northern blots, dot blots			
	The durability and high protein retention of Immobilon PVDF make these membranes the preferred option for sequential detection of multiple targets via antibody stripping and reprobing.							
Selection Criteria	• Target protein expressed over a broad dynamic range • Wide range of molecular weights		• Target protein <20 kDa • Very low abundance target protein • Low affinity antibody		• High-abundance target protein • Wide range of molecular weights		• High-abundance target proteins • Target protein <20 kDa	

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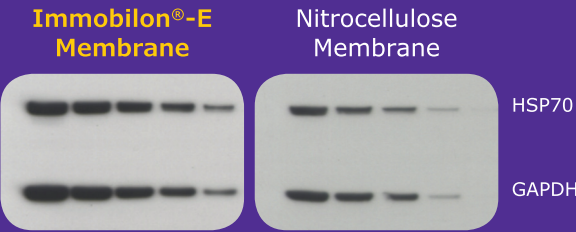
No more Alcohol pre-wet
Alcohol = methanol or ethanol or isopropanol

NEW Immobilon®-E membrane for high sensitivity immunodetection

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The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.
Merck and the Vibrant M are trademarks of Merck KGaA, Darmstadt, Germany. Immobilon is a registered trademark of Merck. Copyright © 2017 Merck KGaA. All Rights Reserved. All Rights Reserved. Lit. No. MK_FL1158EN 2017 - 07460 10/2017

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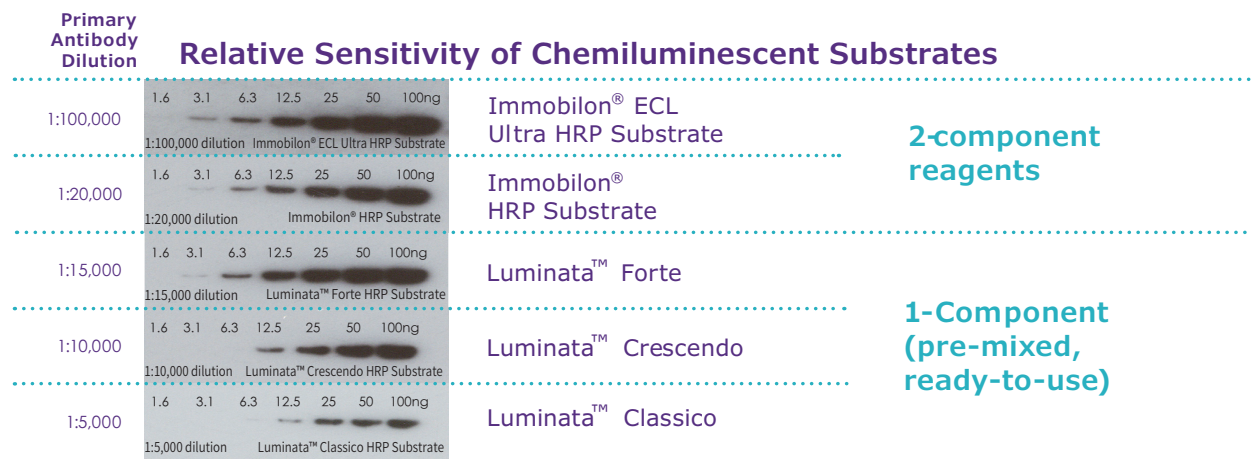
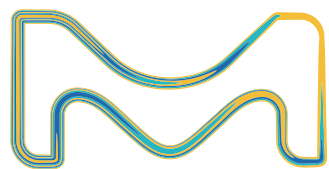
Improve your signal in your next Western blot.
sigma-aldrich.com/westernblot

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and the Vibrant M are trademarks of Merck KGaA, Darmstadt, Germany. Immobilon is a registered trademark of Merck KGaA, Darmstadt, Germany and its affiliates. Copyright © 2016 Merck KGaA. All Rights Reserved.

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ReBlot™ Plus Western blot recycling kit

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What is ReBlot™ Plus?


- Re-Blot™ Plus 溫和型 Stripping Solution - Provides good results on both nitrocellulose and PVDF membranes.
- Re-Blot™ Plus 強力型 Stripping Solution - Performs when membranes with high signal are to be stripped, or use when Re- Blot™ Plus Mild treatment is not sufficient.

Key Benefits


- β-Mercaptoethanol-free to avoid pungent odors
- Room temperature stripping in just 15 minutes
- Fast reuse of blots for multiple antibody probings
- Non-acidic, for reduced risk of protein degradation (such as in Edman degradation)

SNAP i.d.® 2.0 System

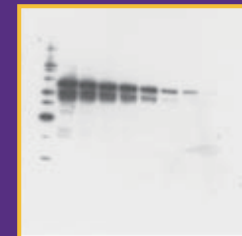
HSP70

Stripped Blot

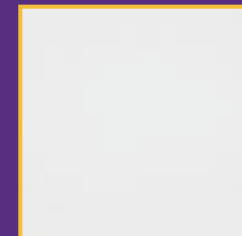GAPDH

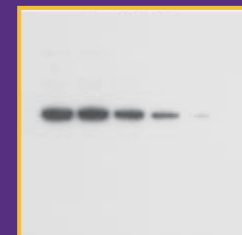ReBlot™ efficiently strips blots to allow for fast reprobing with different antibodies.

Two-fold dilutions of A431 lysate were resolved by SDS-PAGE then transferred onto Immobilon® -P. The blot was initially probed with HSP70 (1:8,000, Cat. No. H5147) (top) using the SNAP i.d.® system. Following stripping using ReBlot™ Plus Strong for 15 minutes (middle), the blot was reprobed with anti-GAPDH (1:8,000, Cat. No. MAB374) (bottom).

ReBlot's™ ability to efficiently strip the blot led to a clean GAPDH blot, even though both primary antibodies share the same anti-mouse secondary antibody.

Immobilon[®] NOW

Measurement
marks on the lid.
No ruler needed.

Immobilon
NOW Dispenser[®]